Miltenyi Biotec

MACSQuant[®] X

User manual

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No part of this user manual may be reproduced, stored in a retrieval system, transmitted, published, or distributed in any form or by any means, electronically, mechanically, by photocopying, microfilming, recording, or otherwise, without the prior written consent of Miltenyi Biotec; however, notwithstanding the foregoing, the owners of the MACSQuant[®] X may make copies solely for purposes of training personnel in the use and servicing of the unit within their business or organization. Maximal care has been taken by Miltenyi Biotec in the preparation of this user manual. However, Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this document. The information in this document is provided "as is" without warranty of any kind, and is subject to change without notice. MACS, the MACS logo, MACSQuant and MACSQuantify[™] are either registered trademarks or trademarks of Miltenyi Biotec GmbH and/or its affiliates in various countries worldwide. All other trademarks mentioned in this document are the property of their respective owners and are used for identification purposes only. Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for therapeutic or diagnostic use.

MACSQuant[®] X

User manual

Version 1

Original instructions

Miltenyi Biotec GmbH
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macs@miltenyibiotec.de
www.miltenyibiotec.com

Read the user manual before using the instrument

Before using the MACSQuant[®] X, read the chapter **Important safety information** and all other information contained in this user manual, including any safety and operating instructions. Pay special attention to all warnings displayed on the instrument. Failure to read and follow these guidelines could lead to improper or incorrect usage and result in damage to the instrument. Improper usage could also cause severe personal injury, death, unpredictable results, instrument malfunction, and premature wear to components shortening the lifetime of the instrument. Such actions may void your warranty. Keep the user manual and any other safety and operating instructions provided with the instrument in a safe place accessible to all users for future reference.

If you have a serious concern regarding the safe use of your instrument, please contact your authorized Miltenyi Biotec service provider or call Miltenyi Biotec Technical Support.

Content

Important safety information	7
Wichtige Sicherheitshinweise	15
Información importante de seguridad	23
Informations de sécurité importantes	31
Importanti istruzioni di sicurezza	39
1 Introduction	47
1.1 Key features of the MACSQuant® X	48
2 Installation	49
2.1 Components included in the delivery	49
2.2 Consumables	50
2.3 Accessories	50
2.4 Service contracts	50
2.5 Unpacking the MACSQuant® X	51
2.6 Assembly	52
2.6.1 Install the fluid bottles	52
2.6.2 Install the Single tube rack	54
2.6.3 Install the MACSQuant X Orbital Shaker	54
2.6.4 Connect the instrument to the power supply	54
2.6.5 Install the webcam	54
3 Operating the instrument	57
3.1 Switching on the instrument	57
3.2 Shutting down	57
3.2.1 Manual shutdown	57
3.2.2 Automatic shutdown	57
3.3 Long-term storage	58
4 Calibration	59
4.1 Calibrating the uptake unit	59

4.1.1	Calibrating the robotic needle arm	59
4.1.2	Calibration of the Single tube rack	61
4.1.3	Calibration of the MACSQuant X 5 Rack	61
4.1.4	Calibration of the 96-well plate	62
4.1.5	Calibration of the 384-well plate	62
4.1.6	Calibration of the MACSPlex Filter Plate	63
4.1.7	Calibration of custom 96 and custom 384 racks	63
4.1.8	Calibration of Universal Reagent Rack	63
4.2	Calibrating the rack detection	64
4.2.1	Single tube rack detection	64
4.3	Volume calibration	64
4.4	Tube length calibration	67
5	Hardware monitor	69
5.1	Hardware monitor	69
5.1.1	Fluidics	69
5.1.2	Sample uptake unit	69
5.1.3	Lasers and detectors	70
6	Live Support	71
7	Maintenance	73
7.1	Decontaminating the instrument	73
7.2	Exchanging fluid bottles	73
7.3	Exchanging hydrophobic air filters	73
7.4	Cleaning the washing station	73
7.5	Cleaning the uptake needle	74
7.6	Exchange the sample needle	74
7.7	Exchange the sample tubing	76
7.7	Exchange the needle sealing	79
8	Troubleshooting	83
8.1	Syringe pump leakage	83

8.2 Air bubbles during measurement / no event acquisition	83
8.3 Excessive debris is present in acquisition	83
8.4 Touchscreen remains dark	83
9 Technical data and specifications	85
9.1 EC Declaration of conformity	88
9.1.1 Additional international conformity	88
10 Technical support	91
11 Limited warranty	93

Important safety information

The instrument is designed for safe use when installed correctly, operated by trained personnel, operated in accordance with general safety practices and the instructions in this user manual.

Warnings and precautions

The guidelines in this section explain the potential risk associated with operating this instrument and provide important supplemental safety information to minimize the risks. Follow the instructions carefully to protect yourself, others, and the equipment from potential hazards and create a safe work environment. Use this instrument only as specified by the manufacturer to avoid damage of equipment and injury to personnel. Always follow local working area safety instructions and laboratory policies as well as standards for health, safety and prevention of accidents. Contact the local authorities governing electrical power supply, building constructions, maintenance, or safety for more information about the safe installation and operation of the instrument.

Hazard levels

Signal words are used to identify safety and property damage messages. The following signal words are used to throughout this user manual.

or **WARNING!** indicates a potentially hazardous situation, which, if not avoided, could result in death or serious injury.

or **CAUTION!** indicates a potentially hazardous situation, which, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices.

Safety symbols

The following symbols are used to highlight conditions that could cause injury to personnel.

Safety alert. Risk of danger. The documentation needs to be consulted in all cases where this safety alert symbol is used on the instrument, in order to find out the nature of the potential hazard and any actions that have to be taken.

Risk of electric shock.

Risk of crushing and shearing of bodily parts.

Hazardous optical radiation.

Hazardous laser radiation.

Biohazard. Risk of contamination if potentially dangerous biological material is used.

 Protective conductor terminal. Symbol is attached on the inside of the instrument. Information for service personnel.

 On (Power supply). Off (Power supply).

Safety labels

Please notice the hazard points and safety symbols of the instrument. All safety labels and safety markings must be kept clean and legible. Inspect the safety labels and safety markings regularly and replace them if not legible or identifiable from a safe viewing distance. Contact Miltenyi Biotec for replacement labels.

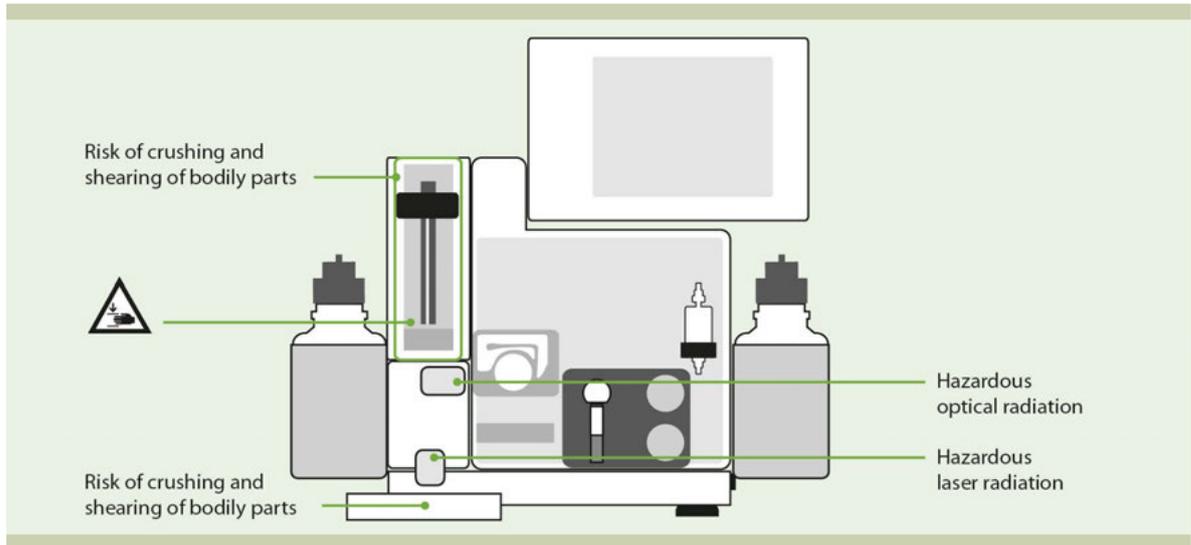
Figure 1: Hazard areas and safety symbols on the front of the MACSQuant® X.

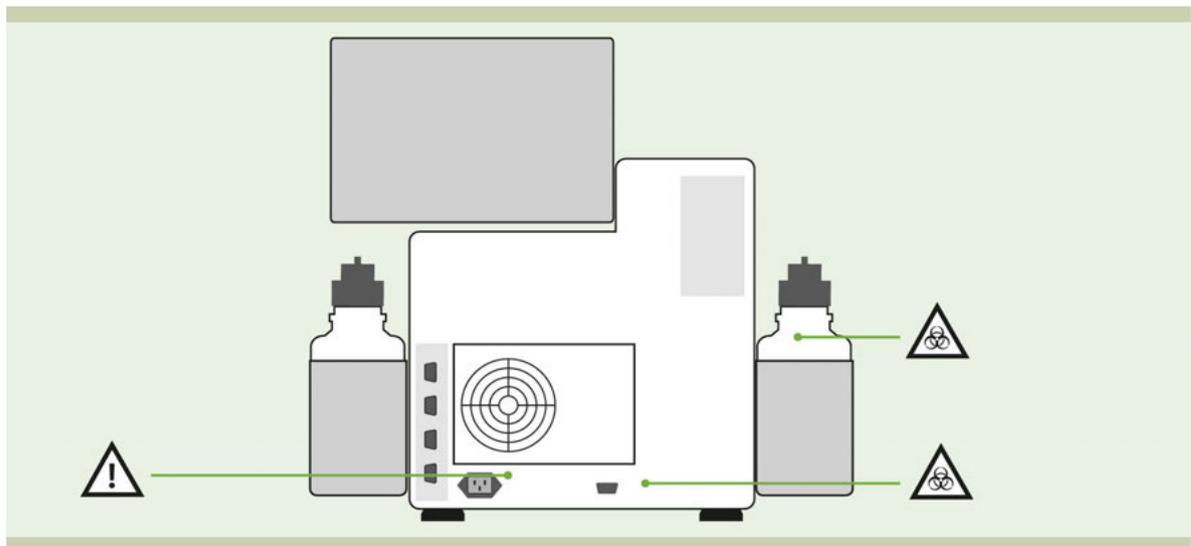
Figure 2: Hazard areas and safety symbols on the rear side of the MACSQuant® X.

General safety instructions

Safe operation of the instrument is no longer ensured if the instrument is not working properly and/or the displayed instructions or messages advise you to contact Technical Support. Immediately switch off and unplug the instrument from the power outlet and contact an authorized Miltenyi Biotec service provider or Miltenyi Biotec Technical Support.

Electricity and fire hazards

Potential risks arising from electrical devices include electric shock, short, and overheating. Electric shock may lead to severe personal injury or death. An electric short and overheating may lead to the spread of fire. Electronic equipment might emit sparks, which could ignite combustible vapors or material resulting in explosion or fire.

Do not use the instrument in areas classified as hazardous locations; for example, in oxygen-laden environments. If flames or smoke appear, immediately switch off the power supply, unplug the instrument from the electrical outlet, and contact an authorized Miltenyi Biotec service provider or Miltenyi Biotec Technical Support.

The instrument housing is designed to lower the risk of electric shock, short, and spread of fire.

Except for the front access covers, do not remove or penetrate any cover, nor any cover of other accessory hardware supplied by Miltenyi Biotec. Only authorized personnel is allowed to remove any other cover of the instrument. Never push an foreign object through an opening into the instrument.

Do not use the instrument if

- it has been opened or disassembled,
- it has been dropped or damaged,
- it has damaged or broken parts,
- it has a damaged power cable,
- an object has been entered through the ventilation slots,
- an object has been dropped into the instrument.

The instrument is intended for indoor use only. If liquids enter the instrument, it may lead to electric shorts, shock, or fire. Take special care while handling liquids.

Do not allow liquids to enter the interior of the instrument. Protect the instrument against accidental spillages and splashes. Clean up spillages immediately. Do not operate the instrument if liquids have entered the instrument. Do not use the instrument in a wet and damp location. Avoid areas with high humidity or condensation. Moving the instrument from a cold environment – such as a cold room at 5°C – to room temperature may cause condensation inside the instrument. In such cases, wait for the instrument to dehumidify before operating it. Before cleaning, unplug the instrument. Do not use liquid or aerosol cleaning agents, always use a damp cloth.

Protect the instrument from overheating. Do not use the instrument in areas with more than 25°C. Ensure adequate air circulation in the room when the instrument is operated. Allow sufficient air circulation around the instrument – at least 15 cm on all sides – during operation. **Do not place the instrument near radiators, heat registers, stoves, or other equipment (including amplifiers) that generate heat. Protect the instrument from direct sunlight. Do not cover the slots and openings of the instrument as these are provided for ventilation.** Do not place the instrument in a built-in rack or similar confined spaces unless the space has been specifically designed to provide proper ventilation.

The instrument is equipped with a three-wire electrical grounding-type plug that has a third pin for grounding. This plug fits in a grounded power outlet only. This is a safety feature.

Do not try to insert the plug into a non-grounded power outlet. If you cannot insert the plug into the outlet, contact your local electrician to replace the outlet.

Only peripheral devices that comply to UL 60950 are allowed to be connected to the RS232 connector labeled COM1.

Only the 2D code reader that was delivered with the instrument should be connected to the RS232/BCR port. External laser devices connected to the RS232/BCR port have to comply with the standard IEC 60825-1.

In order to ensure electromagnetic compliance, connect peripheral devices only with double-shielded cables and only use the power cable and the keypad included in delivery. Use the keypad with MACSQuant X models only.

Ensure that the main switch as well as the plug of the power cable are easily accessible and located as close to the operator of the instrument as possible. If it is necessary to disconnect the power supply, unplug the cable from the power outlet.

The instrument should only be operated from a power source that meets the specifications mentioned on the product's electrical ratings label. If you have questions about the type of power source to use, contact your authorized Miltenyi Biotec service provider or local power company. Do not use extension cords or power strips. Do not overload an electrical outlet.

Mechanical hazards

Moving and revolving parts are potential mechanical hazards.

Do not open the front access covers while the instrument is in operation. Do not obstruct the movement of the robotic needle arm and accessory hardware during operation. There is a risk of puncturing, crushing and shearing bodily parts. Keep fingers etc. away from all moving parts of the instrument or accessory hardware. Do not circumvent any safety measures or devices. Do not touch fluid pumps or adjust the tubing while the instrument is in operation. Always switch off the instrument before adjusting any part of the fluidic system. Always stop or abort a procedure before handling accessory hardware, e.g., MACSQuant X Orbital Shaker, or loading/removing tubes from the tube rack.

Optical radiation hazards

The MACSQuant X is classified as a Class 1 M laser product per standard IEC 60825-1:2014.

The MACSQuant X is equipped with three Continuous-Wave Lasers (laser class 3B). These lasers are secured by a protective housing.

The instrument is equipped with four vertical cavity surface emitting lasers (VCSEL) for automated rack detection (Class 1M). The radiation is not visible.

Invisible laser radiation. Do not expose users of telescopic optics.

The instrument is equipped with a 2D Code Reader and with powerful light emitting diodes (LEDs) for illuminating the reading area. Powerful LEDs are also used to illuminate the fluid bottles as well as the waste bottle. According to the international standard IEC 62471, this lamp system has an exposure hazard value (EHV) of 0.91 and is in excess of the Exempt Risk Group.

Do not look directly at laser or LED radiation or reflected laser or LED radiation from a mirrored surface. Otherwise, eye injury may result. Do not direct the laser beam at others.

Be careful of the path of the light beam or reflection from a mirrored surface. Take care during installation of the instrument that the path of the light beam is not at the same height as the human eye during operation. Do not allow water, oil, dust or other foreign substances to stick to the 2D Code Reader aperture window. This may cause reading errors. Be sure to stop the laser emission before cleaning the scanner. Otherwise, exposure to the light may cause eye injury. Use a soft, dry cloth to wipe any substances from the scanner. Do not use alcohol or other cleaning substances.

Do not disassemble, modify, or remove the installed laser or LED radiation sources. The laser or LED radiation sources do not automatically stop emitting when disassembled. Radiation of disassembled units may lead to eye injuries.

	Output power	Pulse time	Wavelength
MACSQuant X			
Rack Detection	3.3 mW	215 µs	850 nm
Internal Laser 1	30 mW	Continuous	488 nm
Internal Laser 2	40 mW	Continuous	404 nm
Internal Laser 3	20 mW	Continuous	640 nm

Table 2.1: Laser output power, pulse time, and wavelength.

Chemical and biological hazards

Depending on the biological material used, contamination infection may lead to severe personal injury or death. All clinical samples must be considered potentially infectious.

If biohazardous material is or has been used, the operator must choose and wear personal safety equipment as indicated in the warnings and precautions for the particular substance. The above safety precaution also accounts for any hazardous chemical, including toxic or corrosive chemicals, acidic or radioactive substances that may be present in the sample.

Wear protective gloves and clothing, as well as safety glasses to prevent contact of the substance with skin and eyes. Also protect mouth and nose as aerosols might leak from the system. Defective or inadequate safety equipment is hazardous. The instrument must be operated in a safety hood if hazardous or unknown material is being processed. If hazardous material has been used or spilled, care must be taken to thoroughly decontaminate the instrument. It is strictly prohibited to continue to handle contaminated accessories or parts of the instrument.

All liquid and solid waste must be considered hazardous and, must be therefore handled taking universal laboratory precautions. Waste disposal must be in accordance with any local regulations.

Before disposal, decontaminate all tubes and other consumables that were in contact with biohazardous material. Liquid waste must be autoclaved or decontaminated using a disinfectant that is appropriate for the specific pathogen, e.g., 10% bleach, isopropyl alcohol, or 70% ethanol.

Failure of parts containing biohazardous material or liquids that have been in contact with such materials could cause a hazard.

Always inspect the fluidics system (complete tubing set, reservoirs, bottles and their closures, valves, columns, diluters, peristaltic pumps, and uptake needle) before switching on the instrument. If leakage has been detected, replace all damaged parts before switching on the instrument. If damaged part cannot be replaced, unplug and do not use the instrument.

If using reusable steel tips, the waste container may contain potentially infectious material. Label the container properly and dispose of the liquid waste in accordance with your local regulations.

Servicing and transportation

Improper or incorrect servicing or repair of the instrument can cause hazards to users, lead to unpredictable results, cause instrument malfunctions or damage, as well as premature wear and tear and reduced life of the instrument. It may also void your warranty.

Unless otherwise specifically noted in this user manual or other Miltenyi Biotec documentation, do not service the instrument yourself. Servicing and repair must be performed by qualified service personnel.

When replacement or spare parts are required, ensure that the service provider uses only genuine Miltenyi Biotec parts, or third-party parts specified and recommended by Miltenyi Biotec. Using unauthorized parts can cause malfunctions of the instrument and impair results. Miltenyi Biotec does not honor any warranty or accept any responsibility for instrument failure or damages resulting from the use of inappropriate parts. After completing any service or repair work, ensure that your authorized Miltenyi Biotec service provider performs all safety checks as required by the repair procedure to ensure that the instrument is operating correctly.

Only use options and upgrades recommended by Miltenyi Biotec. Inquire with your local Miltenyi Biotec representative about Miltenyi Biotec's extensive instrument service and support arrangements, or refer to www.miltenyibiotec.com/support.

The instrument should be transported with care in packaging specified by Miltenyi Biotec. Internal damage can occur if the instrument is subjected to excessive vibration or is dropped. If the instrument needs to be shipped back to the manufacturer for service, decontaminate the instrument to remove any hazardous material prior to shipment. If you have questions regarding proper decontamination or shipment, contact Miltenyi Biotec Technical Support for assistance.

Disposal

Waste of Electrical and Electronic Equipment (WEEE) customer information

Please dispose of your end-of-life Miltenyi Biotec products in accordance with the applicable WEEE and hazardous waste disposal legislation, which may differ by country or region.

Electrical and electronic equipment may contain hazardous substances which may have a serious detrimental effect on the environment and/or human health. That is why all equipment must be specifically collected and treated by designated waste facility centres and by qualified WEEE compliance schemes. By ensuring that you dispose of your unwanted electrical and electronic equipment according to the applicable WEEE and hazardous waste disposal legislation, you are helping to preserve our natural resources and protect human health.

Miltenyi Biotec is committed to protecting the environment. Miltenyi Biotec offers product end-of-life return programs in many countries, and partners with licensed WEEE compliance schemes throughout the world. Miltenyi Biotec lets you recycle your end-of-life Miltenyi Biotec equipment free of charge. The terms and availability of this offer vary by geography because of differences in regulatory requirements. Please note that, depending on the type and use of your equipment, additional requirements may apply.

For more information, or if you wish to dispose of your end-of- life Miltenyi Biotec equipment, please contact your local Miltenyi Biotec representative or Miltenyi Biotec Technical Support.

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1

Introduction

MACSQuant® Instruments are compact flow cytometers that combine multisample and multiparameter sample analysis with unrivaled ease of use. With processing rates of up to 15,000 events per second, they facilitate also absolute quantitation of cell populations. Miltenyi Biotec is leading provider of comprehensive solutions for flow cytometry.

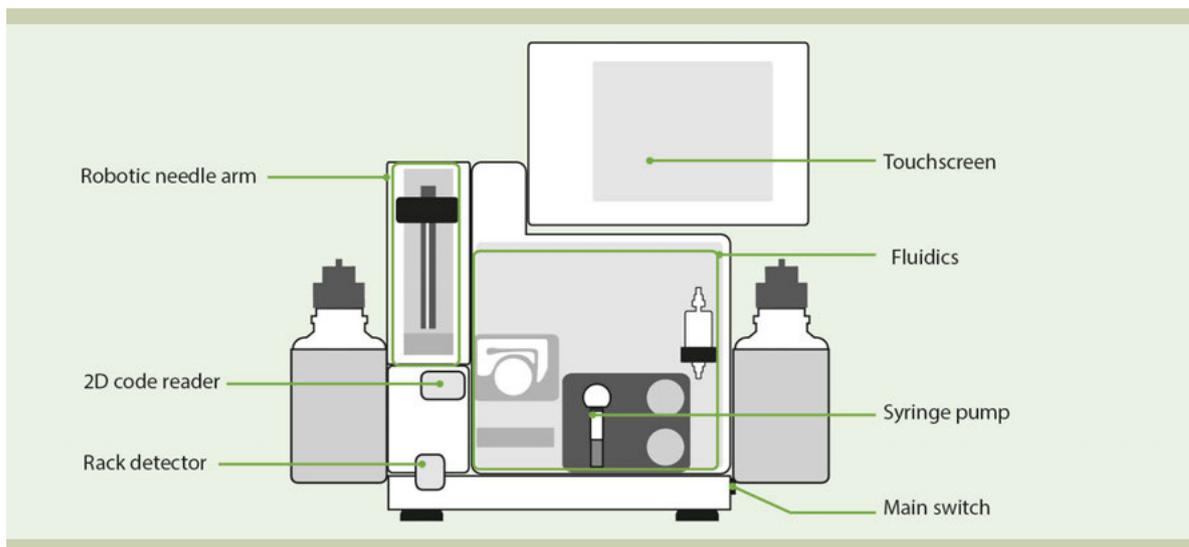
Figure 1.1: Front view of the MACSQuant® X.

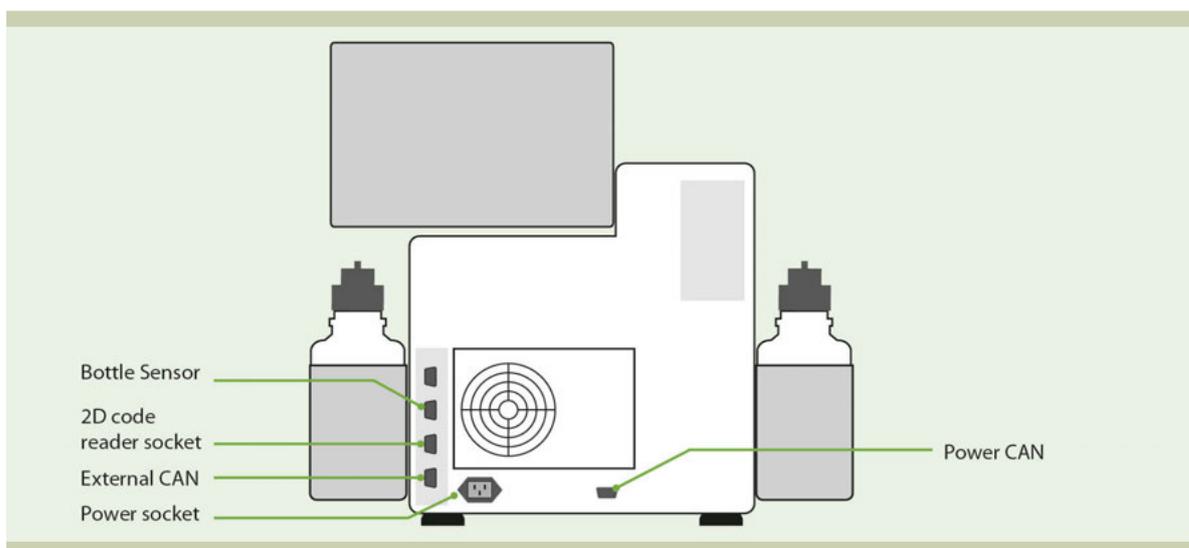
Figure 1.2: Rear view of the MACSQuant® X.

1.1 Key features of the MACSQuant® X

The MACSQuant X has been engineered from the ground up as a compact flow cytometer made to provide maximum speed and reliability in high throughput screening. It allows you to get more data, in less time, with less effort. Perfect for laboratories that want to optimize time while enjoying the certainty of getting reliable results every single run, the MACSQuant X has been designed to meet high-throughput cell analysis needs for the most challenging large-scale screening settings.

High-throughput

Process and analyze up to 10 parameters in 384- and 96-well formats.

Speed

Run an entire 96-well plate in only 15 minutes and a whole 384-well plate in under 60 minutes.

Reliability

Get consistent results every experiment regardless of sample type and volume.

Flexibility

Easily and immediately switch between tubes and plates at the click of a button.

Automation

Take advantage of a truly hands-free instrument, which can be easily integrated into liquid handlers.

2

Installation

If there is any damage, do not use the instrument but contact your local Miltenyi Biotec representative or Miltenyi Biotec Technical Support. The operation of a damaged instrument may lead to the exposure of mechanical hazards, optical radiation hazards, electric shock, or the spread of fire. Read the chapter **Important safety information** before continuing with installation and assembly. Read through the following instructions carefully before commencing the installation procedure.

Before opening the transportation box, check for any visible external damage to the box. Check also to see if the shock and position indicators (if present) suggest incorrect transportation of the instrument.

The MACSQuant[®] X is designed to fit perfectly onto benchtops, into standard size laminar flows, into safety cabinets, or in connection to fully automated liquid handling systems. It should be installed on a stable, flat and vibration-free surface. The operating environment should be dust-free, sufficiently ventilated, and free from sources of electromagnetic radiation. In order to ensure a flat surface in laminar flow hoods, the MACSQuant[®] X can be placed on a MACS Laminar Hood Plate.

2.1 Components included in the delivery

The following components are included in the delivery of the X (# 130-105-100):

Component	Position in transportation box
MACSQuant [®] X	Main storage compartment
Fluid sensor cable module	Top storage compartment
Power cable	Top storage compartment
Fluidics tubing (6x)	Top storage compartment
Bottle closure with sensors (4x)	Top storage compartment
Single tube rack	Top storage compartment
User manual	Top storage compartment
Empty bottles with closures	In bottle holders
Keyboard	Front compartment
Webcam	Top storage compartment
Hydrophobic Air Filters	Top storage compartment
MACSQuant Calibration Beads	Top storage compartment
MACSQuant Starting Buffer Kit	Extra packaging

2.2 Consumables

Component	Description	Order no.
Hydrophobic Air Filter	To vent fluid bottles	130-090-385
MACSQuant Calibration Beads	For instrument calibration	130-093-607
MACSQuant Running Buffer	6 × 1.5 L	130-092-747
MACSQuant 16× Running Buffer	1 × 1.25 L	130-111-562
MACSQuant 16× Running Buffer	6 × 1.25 L	130-111-747
MACSQuant Washing Solution	6 × 1.5 L	130-092-749
MACSQuant Storage Solution	6 × 1.5 L	130-092-748
MACSQuant Starting Buffer Kit	4 × 1.5 L Running Buffer 1 × 1.5 L Washing Solution 1 × 1.5 L Storage Solution	130-094-190
MACSQuant Washing & Storage Solution Kit	3 × 1.5 L Washing Solution 3 × 1.5 L Storage Solution	130-092-801
MACS Bleach Solution	1 × 1.0 L	130-093-663
MACS Comp Bead Kit, anti-human Igk	for 100 tests	130-104-187
MACS Comp Bead Kit, anti-mouse Igk	for 100 tests	130-097-900
MACS Comp Bead Kit, anti-rat Igk	for 100 tests	130-107-755
MACS Comp Bead Kit, anti-REA	for 100 tests	130-104-693

2.3 Accessories

Complementary components can be ordered separately.

Component	Description	Order no.
MACSQuant X Orbital Shaker	Use the MACSQuant X Orbital Shaker with the MACS Universal Reagent Rack and the MACSQuant X 5 Rack for 5 mL conical tubes.	130-112-558
Buffer Supply Station	To connect larger fluid containers	130-101-841
Racks	Universal Reagent Rack MACSQuant X 5 Rack	130-115-722 130-112-413


2.4 Service contracts

For information about available service contracts for the MACSQuant X, please visit www.miltenyibiotec.com/.


2.5 Unpacking the MACSQuant[®] X

The cardboard packaging constitutes of three pieces that are held together by tensioning straps.

- 1 Remove the tensioning straps.
- 2 Lift the top cardboard cover. **Note:** Do not use the handles on the middle cardboard piece to open the packaging at this time.
- 3 The top storage compartment holds the MACSQuant[®] X user manual, cardboard boxes and bags, which contain components included in the delivery as well as any complementary components.
- 4 Take out all boxes and bags from the top storage compartments and remove the foam piece.
- 5 Holding its handles, lift and remove the middle cardboard piece. Stow it aside. **Important:** Do not lift the foam upwards as this might damage the instrument. The foam can be folded back for removal.
- 6 Carefully fold back foam piece on the side where the needle arm of the instrument is located. Be sure not to damage the CF card eject button of the PC (on rear side of the instrument).
- 7 Carefully slide the protection foam off the instrument in the opposite direction of the needle arm and along the instrument display.
- 8 Lift the MACSQuant[®] X by holding it at the base of the orange bottle holders at both sides of the device. **CAUTION!** Two persons are required to lift the MACSQuant[®] X. Take care to avoid personal strain or injury. The instrument is heavier at the front and should be stabilized while lifting it. Miltenyi Biotec accepts no liability for potential injuries sustained during lifting and/or movement of the device.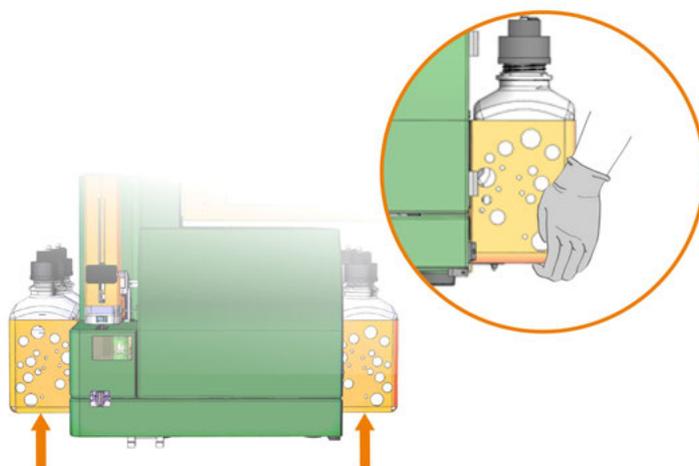

- 9 Transfer the instrument onto a stable working platform and remove the plastic bag surrounding the device. **WARNING!** The instrument requires adequate air circulation for heat exchange and cooling. Read the chapter **Important safety information** to avoid the risk of overheating. Do not place the instrument on an unstable table, cart, stand, tripod, or bracket. As a consequence, the instrument might fall down, which may cause serious bodily harm and/or serious damage to the instrument. Use only a table, cart, stand, tripod, or bracket that can easily support a weight of >50 kg.
- 10 Carefully remove the uptake needle from the foam packaging.
- 11 Place the uptake needle into its guiding in the robotic needle arm.

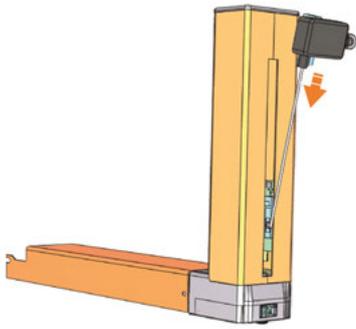
- 12 Remove the protection foam from the waste and storage bottles.
- 13 Remove the stylus from the packaging and place it in the holder attached to the touchscreen.
- 14 Adjust the angle of the touchscreen in an upright position.

2.6 Assembly


2.6.1 Install the fluid bottles

Operating the MACSQuant[®] X requires three fluid bottles for Running buffer, Washing Solution, and Storage Solution. One bottle is required for waste collection. Always operate the instrument with ready-to-use MACS Buffers and solutions only. The bottles, closures, tubing and fluid sensor cables are color-coded for easier handling.


- 1 Install only one fluid bottle at a time. Place the new bottle into the orange bottle holder. Please note the corresponding color coding.
- 2 Unscrew the lid of the fluid bottle and replace it with the appropriate bottle closure delivered with the MACSQuant[®] X.
- 3 Note the color coding of the tubes and connect them to the respective bottle closure. The Running Buffer and the Waste bottle lids have each two connectors for tubes (male and female connector), Washing Solution and Storage Solution bottles have only one.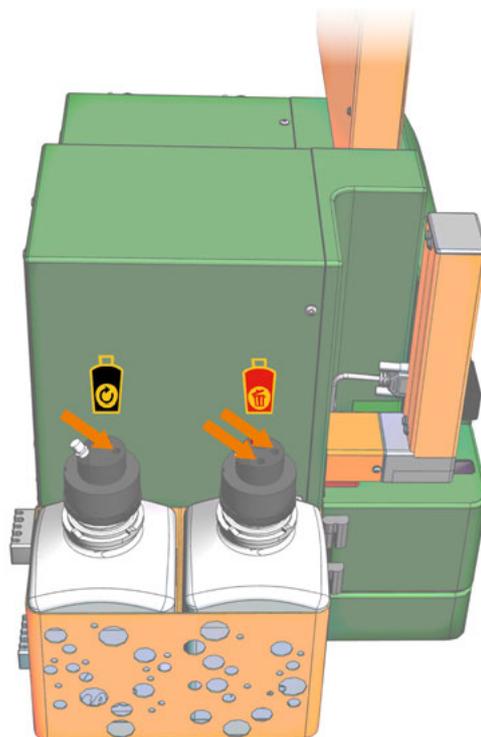

- 4 Connect the 4-end fluid sensor cable to the port labeled **Bottle Sensor** at the back of the MACSQuant® X (**Figure 1.2**) and fasten securely.
- 5 Note the color coding and connect each fluid sensor cable to the respective bottle closure.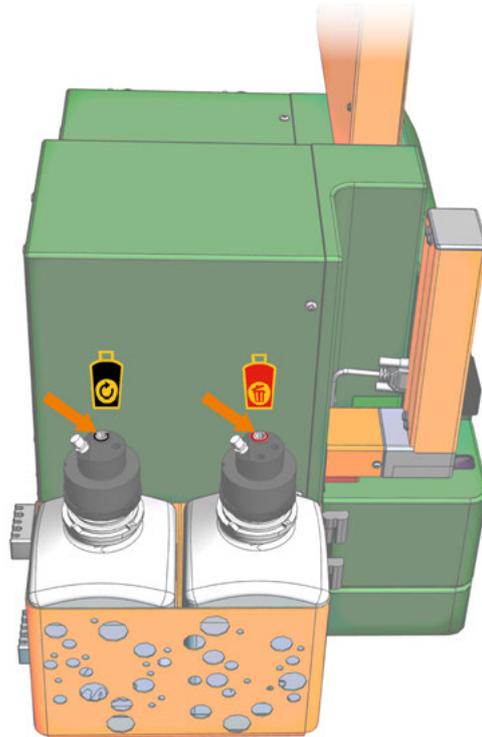

- 6 Connect a Hydrophobic Air Filter to the appropriate connectors on the bottle closures to keep buffers sterile. Avoid any contact with liquids, as this may cause clogging of the filter.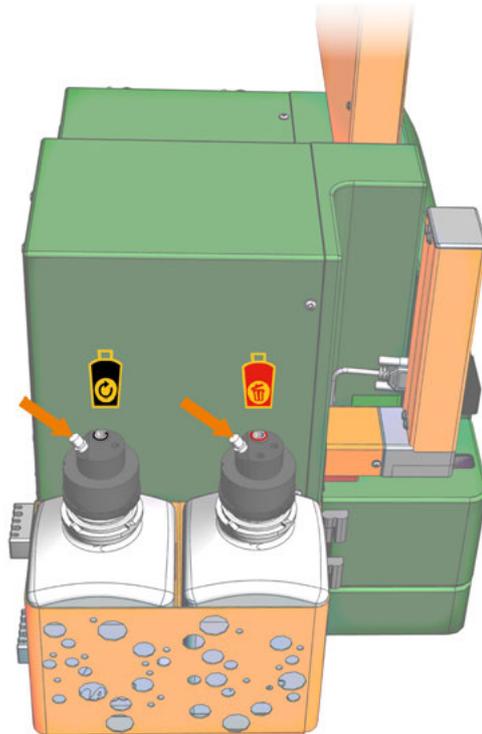


- 7 Remove the protective cap and connect a Hydrophobic Air Filter to the appropriate connector of the fluidics system on the front of the MACSQuant® X.

2.6.2 Install the Single tube rack

The MACSQuant[®] X is delivered with the Single tube rack. The Single tube rack is compatible with all standard flow cytometry tubes (1.5 mL, 2 mL, and 5 mL).

- 1 Gently insert the Single tube rack into the corresponding slots located at the front of the MACSQuant[®] X. The rack should click into place. To remove the Single tube rack, hold it between two fingers and pull gently.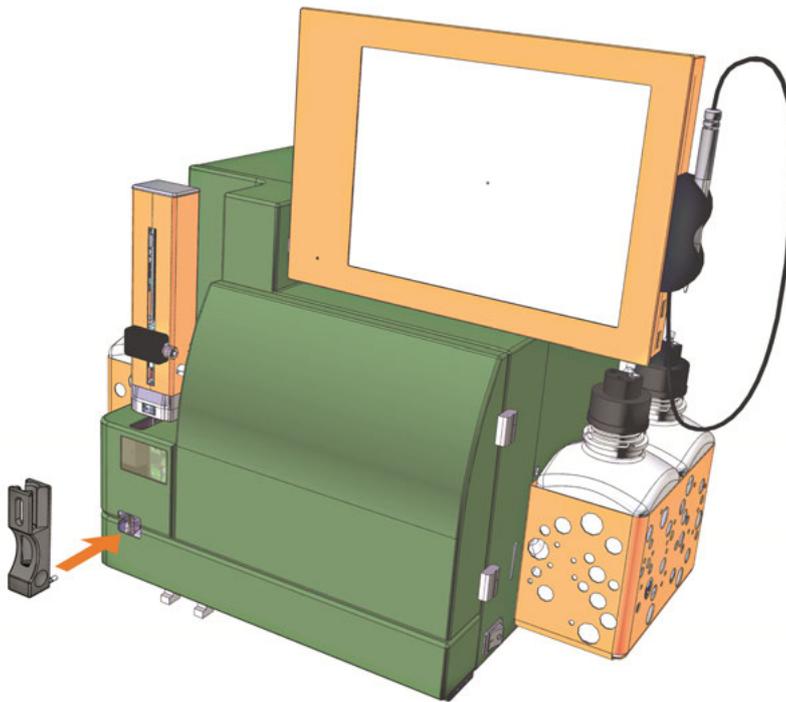

2.6.3 Install the MACSQuant X Orbital Shaker

Refer to the MACSQuant X Orbital Shaker user manual for details on its installation.

2.6.4 Connect the instrument to the power supply

- 1 Connect the power cord to the power socket at the rear side of the MACSQuant[®] X. The main power switch is located on the right side of the instrument (**Figure 1.2**). Ensure that the main switch is in position O (off) before connecting the power cord.
- 2 Attach the cable coming from the 2D code reader to the port labeled **RS232/BCR** at the back of the MACSQuant[®] X and fasten securely.
- 3 Plug in the power cord to a grounded power outlet.
- 4 Switch on the MACSQuant[®] X to turn it into stand-by mode. The MACSQuantify™ Software will automatically power up if you touch the touchscreen.
- 5 Wait until the MACSQuantify™ Software is loaded completely. Connect the keyboard included in the delivery. Connect a mouse if desired.

2.6.5 Install the webcam

- 1 Attach the magnetic webcam holder to the side of the instrument.
- 2 Plug in the webcam cable into a free USB port at the rear side of the instrument.

- 3 Place the webcam into its holder.
- 4 The webcam is recognized automatically.

Note: Unplug the webcam during normal operation of the instrument. Use it only during a Live support session.

3

Operating the instrument

3.1 Switching on the instrument

- 1 Switch on the instrument by using the main switch on the right side of the MACSQuant[®] X. The instrument will be in stand-by mode.
- 2 Touch the touchscreen to power up the instrument.
- 3 For a detailed instruction of how to use the MACSQuantify[™] Software, refer to the MACSQuantify Software User Manual.

3.2 Shutting down

- 1 Shut down the software before switching off the instrument. Shut down the software manually, or enable an automatic shutdown function (refer to section **Automatic shutdown**, page 57).
- 2 Switch off the instrument using the main power switch at the right side of the instrument.

3.2.1 Manual shutdown

- 1 Click the **Main instrument control button** at the upper right corner of the screen.
- 2 Select **Instrument off**. This automatically starts a wash program that lasts for approximately seven minutes. The fluidics system is thoroughly washed with Washing Buffer and finally filled with storage solution.
- 3 At the end of the washing program, the instrument is in standby mode. It can be switched off, or rebooted by touching the touchscreen.

Note: With the **main instrument control button**, the instrument can also be turned into data analysis mode. This will start the washing procedure as well, but allows data analysis in the meantime.

3.2.2 Automatic shutdown

By default, automatic shutdown is enabled. If the instrument is idle for a certain amount of time, it will automatically shut down or, alternatively, switch into **Analysis mode** to elongate diode lifespan of the lasers.

- 1 To modify the automatic shutdown settings, log in as an administrator.
- 2 Go to **Edit > Options (default) > Software > Timers**.
- 3 Set the following parameter:
 - Standby timer, the amount of idle time before shutdown begins.
 - Shutdown timer, the incubation time during the shutdown procedure (5 minutes is recommended)
 - Needle priming time, to avoid that the needle runs dry (20 seconds are recommended)
- 4 Set the **Shutdown default behavior** to either **Analysis mode** or **Instrument off**.

3.3 Long-term storage

- 1 If the instrument is not in use for a longer time, flush the fluidics system every two weeks to prevent blockages.
- 2 Switch on the instrument into acquisition mode. This will initiate the priming of the instrument.
- 3 After the system has been primed and all system checks have been performed, shut down the instrument.

4

Calibration

WARNING Read the chapter **Important Safety Information** in the MACSQuant® X user manual before operating the instrument. When processing infectious, radioactive, poisonous, or any other hazardous liquids, always abide by the necessary safety precautions.


4.1 Calibrating the uptake unit

Before using the MACSQuant® Instrument for the first time, calibrate the hardware. Calibrate also the photomultiplier tubes (PMTs) of the instrument. In order to calibrate the MACSQuant Instrument, the user must be familiar with the MACSQuantify Software. It is highly recommended that hardware calibration is only performed by an administrator who has been trained by Miltenyi Biotec. Refer to the MACSQuant Instrument software guide and/or the help function of the MACSQuant Instrument for an introduction to the MACSQuant Instrument user interface and a detailed description of the instrument settings.


4.1.1 Calibrating the robotic needle arm

During calibration, the robotic needle arm moves between the samples and the dropping station. Ensure that the robotic needle arm can move freely and that no object is obstructing it.


- 1 Click on the **Tools** tab on the sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. A red closed circle indicates that the respective uptake component is not calibrated.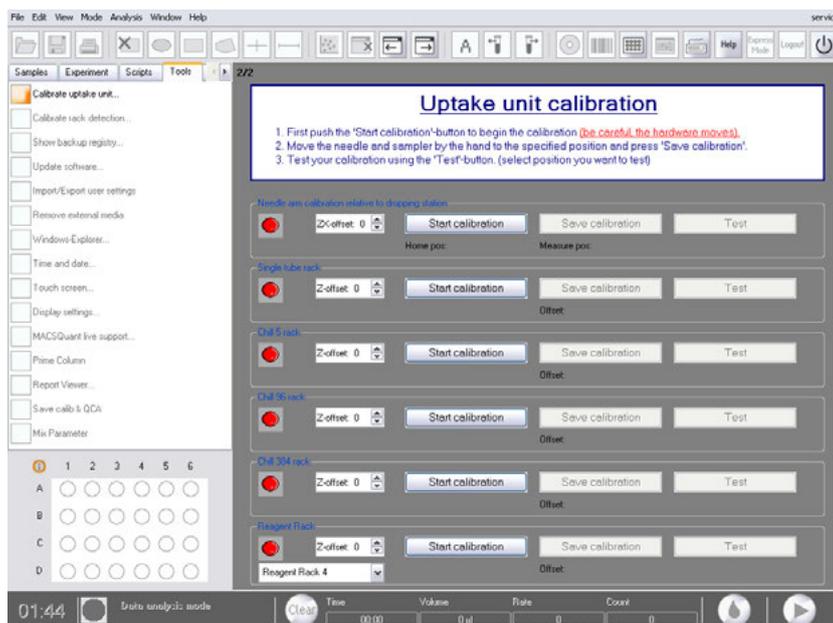

- 2 Click **Start** calibration under the heading **Needle arm calibration relative to dropping station**. The robotic needle arm will move toward the dropping station before being inserted into the port.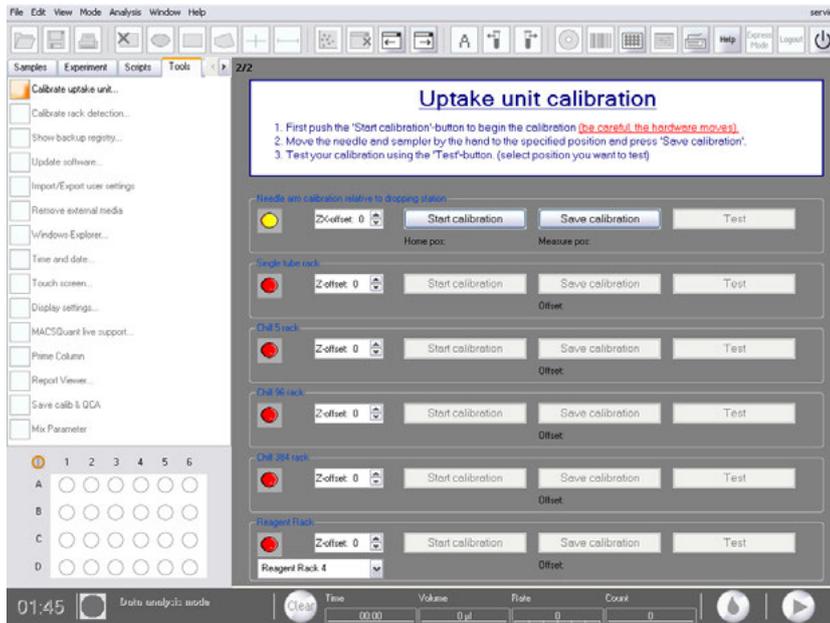

- 3 Manually adjust the robotic needle arm in order to position the uptake needle over the center of the port.
- 4 Gently lower the robotic needle arm into the port until the uptake needle touches the bottom of the port. Do not bend the needle or insert it diagonally.
- 5 Click **Save** calibration.
- 6 The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved.

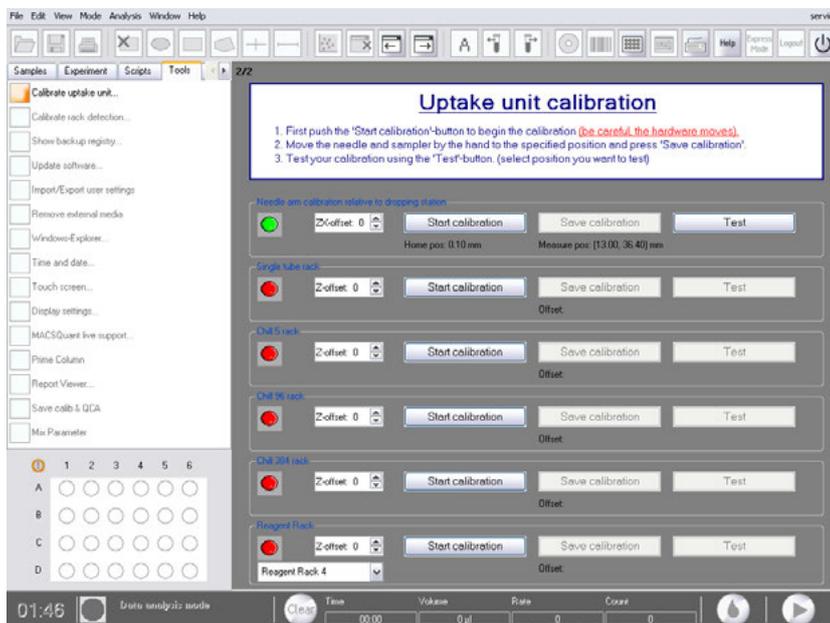4.1.2 Calibration of the Single tube rack

- 1 Attach the Single tube rack with desired sample tube to the instrument as described in **Install the Single tube rack on page 54**.
- 2 Click **Start calibration**. The robotic needle arm will automatically move forward and insert the uptake needle into the single tube. The uptake needle should be positioned at the center of the tube (i.e. equidistant from the tube edges) almost touching the tube bottom (i.e. only a fraction of a millimeter from the bottom of the tube). To check the uptake needle position, gently wiggle the tube to ensure that there is a small amount of movement.
- 3 If this is not the case, carefully adjust the robotic needle arm accordingly.
- 4 Click **Save calibration**. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved.

4.1.3 Calibration of the MACSQuant X 5 Rack

For the correct usage of the MACS Orbital Shaker, please calibrate the MACSQuant X 5 Rack.

- 1 Attach the MACSQuant Orbital Shaker to the MACSQuant Instrument. Refer to the MACSQuant Orbital Shaker user manual for details.
- 2 Click on the **Tools tab** on the sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 3 Click the **Experiment tab** on the sidebar of the MACSQuantify Software and select **Chill 5 rack** from the Rack pull-down menu. If the MACSQuant X 5 Rack is not selected, the MACSQuant Instrument will display a prompt screen.
- 4 Place the MACSQuant X 5 Rack onto the Orbital Shaker and place an empty 5 mL tube in position D6.
- 5 Click **Start calibration**. The robotic needle arm will automatically move forward and insert the uptake needle into the tube in the position D6 of the rack. The uptake needle should be positioned at the center of the tube (i.e. equidistant from the tube edges) almost touching the tube bottom (i.e. only a fraction of a millimeter from the bottom of the tube). To check the uptake needle position, gently wiggle the tube to ensure that there is a small amount of movement.
- 6 If this is not the case, carefully adjust the robotic needle arm accordingly.
- 7 Click **Save calibration**. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved. The MACSQuant Instrument will automatically test selected positions.

4.1.4 Calibration of the 96-well plate


- 1 Click on the **Tools tab** on the Sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 2 Click the **Experiment tab** in the Sidebar of the MACSQuantify™ Software and select the 96-well format from the Rack pull-down menu. If the 96-well is not selected, the MACSQuant Instrument will display a prompt screen.
- 3 Load an empty 96-well plate on the MACSQuant Orbital Shaker.
- 4 Click **Start calibration**. The robotic needle arm will automatically move forward and insert the uptake needle into the well position H12 of the 96-well plate.
- 5 The uptake needle should be positioned at the center of the well (i.e. equidistant from the well edges) almost touching the well bottom (i.e. only a fraction of a millimeter from the bottom of the well). To check the uptake needle position, gently wiggle the plate to ensure that there is a small amount of movement.
- 6 If this is not the case, carefully adjust the robotic needle arm accordingly.
 - a) If the needle is on the left corner of the well, first push the MACSQuant Orbital Shaker from the right, such that the needle is on the right corner of the well. Then proceed with b)
 - b) If the needle is on the right corner of the well, push the MACSQuant Orbital Shaker from the left to reach the center position of the needle in the well. The last movement of the MACSQuant Orbital Shaker must be from the left.
 - c) If the needle is on the rear corner of the well, first move the needle arm to the front, such that the needle is on the front corner of the well. Then proceed with d)
 - d) If the needle is on the front corner of the well, move the needle arm to the front to reach the center position of the needle in the well. The last movement of the needle arm must be from the front.
 - e) Push the needle down, almost touching the well bottom (i.e. only a fraction of a millimeter from the bottom of the well). To check the uptake needle height, gently wiggle the plate to ensure that there is a small amount of movement.
- 7 Click **Save calibration**. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved. The instrument will automatically test selected positions.

4.1.5 Calibration of the 384-well plate

- 1 Click on the **Tools tab** on the sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 2 Click the **Experiment tab** on the sidebar of the MACSQuantify™ Software and select the 384-well format from the Rack pull-down menu. If the 384-well is not selected, the MACSQuant Instrument will display a prompt screen.
- 3 Load an empty 384-well plate onto the MACSQuant Orbital Shaker.
- 4 Click **Start calibration**. The robotic needle arm will automatically move forward and insert the uptake needle into the well position H24 of the 384-well plate. The uptake needle should be positioned at the center of the well (i.e. equidistant from the well edges) almost touching the well bottom (i.e. only a fraction of a millimeter from the bottom of the well). To check the uptake needle position, gently wiggle the plate to ensure that there is a small amount of movement.
- 5 If this is not the case, carefully adjust the robotic needle arm accordingly (refer to section **Calibration of the 96-well plate**, page 62).
- 6 Click **Save calibration**. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved. The instrument will automatically test selected positions.

4.1.6 Calibration of the MACSPlex Filter Plate

- 1 Click on the **Tools tab** on the Sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 2 Click the **Experiment tab** on the sidebar of the MACSQuantify Software and select **MACSPlex Filter Plate** from the **Rack** pull-down menu. If the MACSPlex Filter Plate is not selected, the MACSQuant Instrument will display a prompt screen.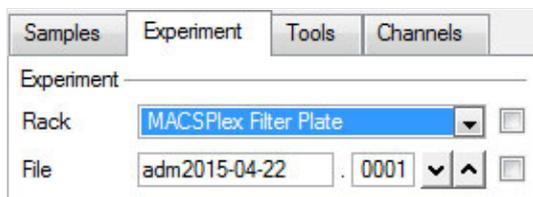

- 3 Place a Chill 96 Rack onto the MACSQuant Orbital Shaker and load an empty MACSPlex Filter Plate.
- 4 Click **Start** calibration. The robotic needle arm will automatically move forward and insert the uptake needle into the rack position H12 of the MACSPlex Filter Plate. The uptake needle should be positioned at the center of the well (i.e. equidistant from the well edges) almost touching the well bottom (i.e. only a fraction of a millimeter from the bottom of the well). To check the uptake needle position, gently wiggle the plate to ensure that there is a small amount of movement.
- 5 If this is not the case, carefully adjust the robotic needle arm accordingly.
- 6 Click **Save** calibration. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved. The MACSQuant X will automatically test selected positions.

4.1.7 Calibration of custom 96 and custom 384 racks

If 96 and 384 plates of different format (e.g. deep well or round bottom) are used, they might have a different distance from the bottom of the MACSQuant Orbital Shaker to the bottom of the well. In order to pick the samples correctly, individual "custom" racks can be defined for each required rack format.

- 1 Click on the **Tools tab** on the sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 2 Click the **Experiment tab** on the sidebar of the MACSQuantify Software and select **Chill 5 rack** from the Rack pull-down menu. If the MACSQuant X 5 Rack is not selected, the MACSQuant Instrument will display a prompt screen.
- 3 Proceed as described for standard 96 or 384 plates (see **Calibration of the 96-well plate on the previous page** or **Calibration of the 384-well plate on the previous page**, respectively).

4.1.8 Calibration of Universal Reagent Rack


- 1 Click on the **Tools tab** on the sidebar of the MACSQuantify™ Software and check the **Calibrate uptake unit** box. Red closed circles indicate that the uptake components are not calibrated.
- 2 Place the Universal Reagent Rack onto the MACSQuant Orbital Shaker.
- 3 Load an empty reagent vial in position 4 (front-most position) of the Universal Reagent Rack.
- 4 Click **Start calibration**. A dialog box will prompt you to ensure that the Universal Reagent Rack is installed correctly and that the Single tube rack is removed.
- 5 Confirm the dialog box by clicking **OK**. The robotic needle arm will automatically move forward and insert the uptake needle into the reagent vial position 4. If this is not the case, carefully adjust the robotic needle arm accordingly.

- 6 Click **Save calibration**. The closed green circle indicates that the calibration is completed. Click **Test** to confirm that the correct coordinates are saved. The MACSQuant Instrument will automatically test selected vial positions.

4.2 Calibrating the rack detection

4.2.1 Single tube rack detection


- 1 Click on the **Tools** tab on the sidebar of the MACSQuantify Software and check the **Calibrate rack detection** box. Red closed circles indicate that the components are not calibrated.
- 2 Click **Start** calibration under the heading **Single tube rack detection**.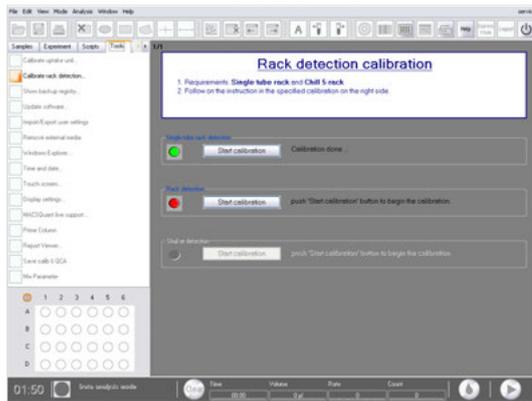


4.3 Volume calibration

After exchange of the sample tubing, the system must be calibrated to this specific tubing in order to correctly set the tube-volume and flow rate. This process is automated as an express mode **Volume Calibration**. Repeat once a month.


- 1 Install either the single tube rack or the MACSQuant X 5 Rack. If using the MACSQuant X 5 Rack, select a single position.
- 2 Select the **Express Mode** of type **Setup** and mode **VolumeCalibration**.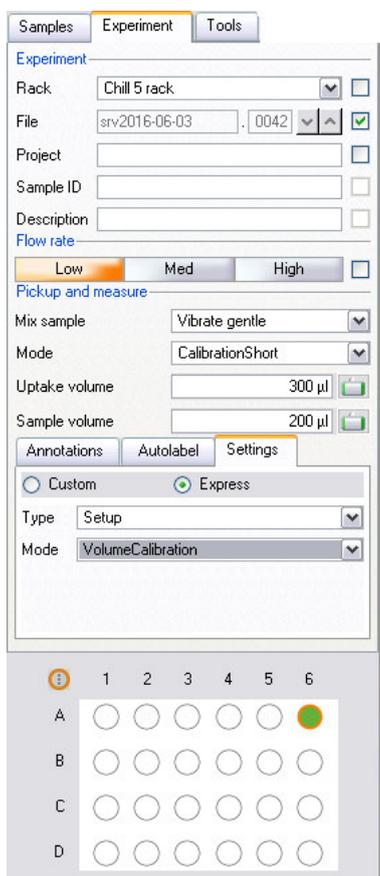

- 3 Click the **Start** button.

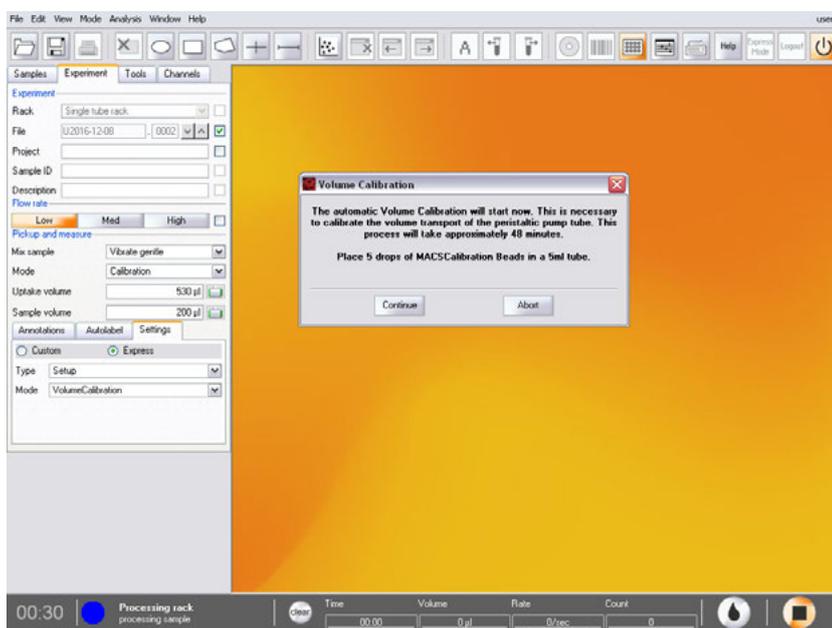- 4 Ensure that buffer bottle is full and the waste bottle is empty. Follow the instruction and click **Continue**.
- 5 After the volume calibration has completed, the status of the calibration is indicated in the upper right corner. **Passed** indicates that the calibration was successful. Otherwise **Failed** indicates that the new tubing could not be identified correctly. In this case, repeat the calibration. If problem persists, contact technical support.

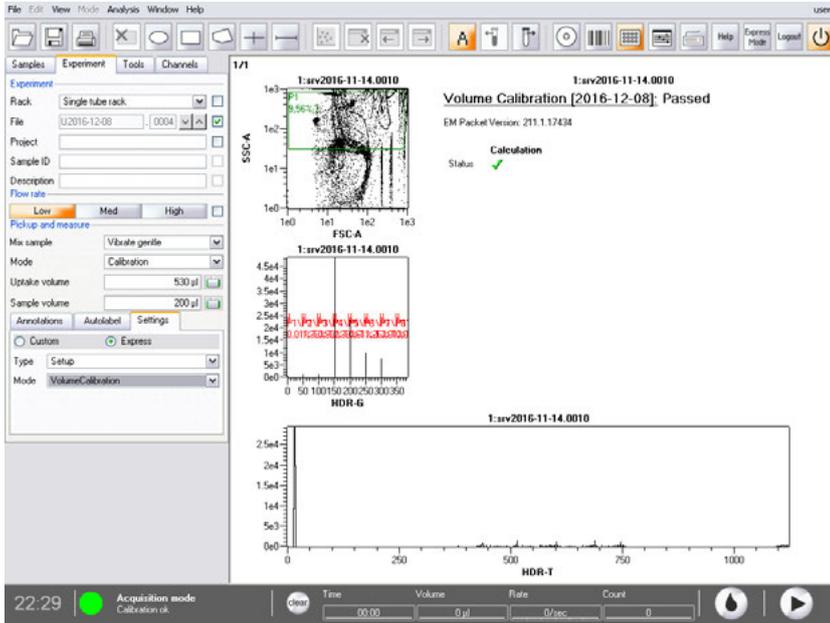4.4 Tube length calibration


The sample tubing is subject to wear, which could affect the tube-volume. In order to renew the teaching process, the Express Mode **Tubing Volume** should be run once a week and will take ~12min.

The Express mode **TubeLengthCalibration** can be performed on the single tube holder or on a single position on the MACSQuant X 5 Rack.


- 1 Select Express-Type **Setup** and Express-Mode **TubeLengthCalibration**.

The screenshot shows the software interface for the MACSQuant X 5 Rack. The 'Experiment' tab is active, showing settings for 'Chill 5 rack' and 'File: srv2016-06-03 . 0042'. The 'Flow rate' is set to 'Low'. Under 'Pickup and measure', 'Mix sample' is 'Vibrate gentle', 'Mode' is 'CalibrationShort', 'Uptake volume' is '300 µl', and 'Sample volume' is '200 µl'. The 'Settings' tab is selected, showing 'Custom' and 'Express' options. 'Type' is 'Setup' and 'Mode' is 'TubeLengthCalibration'. At the bottom, a 4x6 grid of tube positions (A-D, 1-6) is shown, with the tube at position A6 highlighted in green.

- 2 Click the **Start** button.
- 3 Follow the instruction and confirm with **Continue**.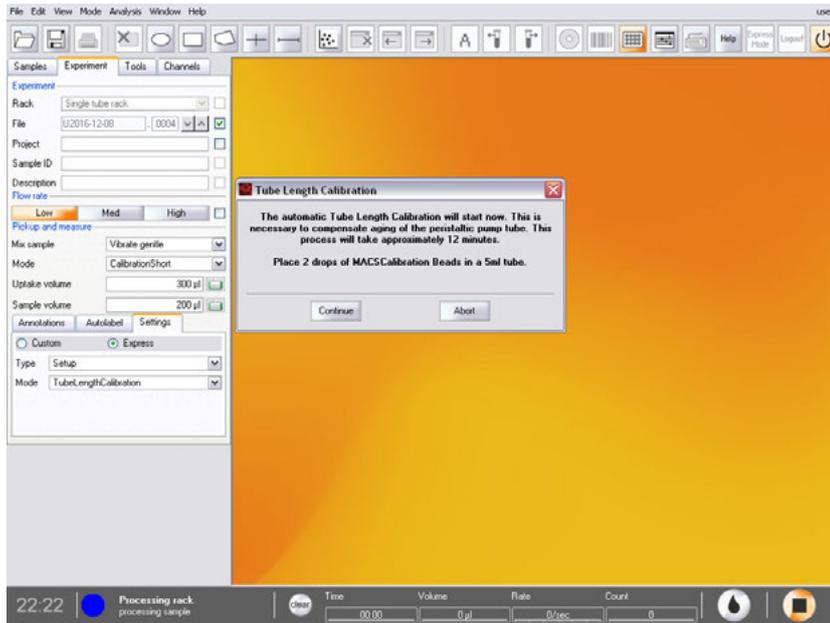

- After the tube length calibration has completed, the status of the calibration is indicated in the upper right corner. **Passed** indicates that the calibration was successful. Otherwise **Failed** indicates that the new tubing could not be identified correctly. In this case, repeat the calibration. If the problem persists, contact technical support.

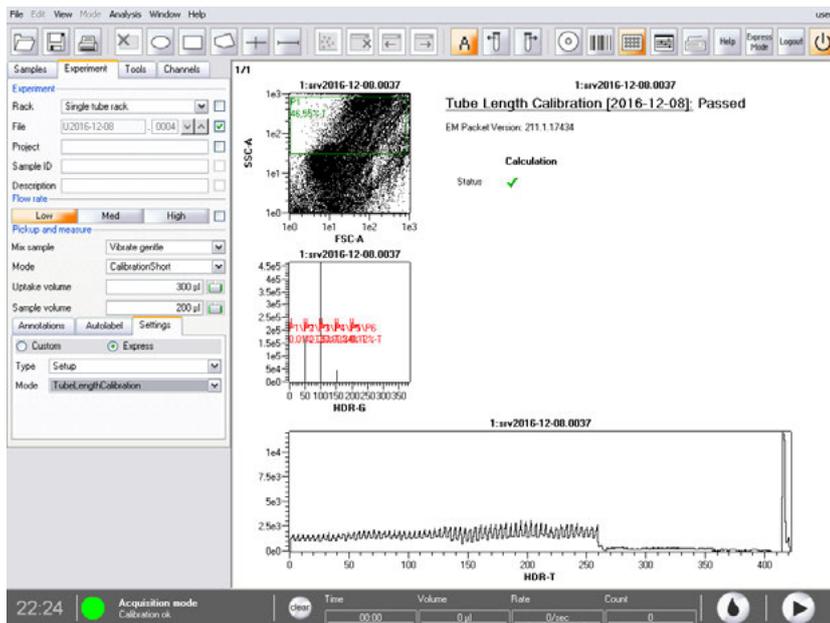5

Hardware monitor

Read the chapter **Important Safety Information** in the MACSQuant® X user manual before operating the instrument. When processing infectious, radioactive, poisonous, or any other hazardous liquids, always abide by the necessary safety precautions.

5.1 Hardware monitor

The hardware monitor can be accessed under the **View menu**. With the hardware monitor the real time status of the MACSQuant Instrument can be assessed, for example, to provide additional information in the case of error messages appearing on the screen. Here you find also the serial number of the MACSQuant Instrument and informations on the photomultiplier tubes (PMT).

5.1.1 Fluidics

The fluidics tab in the hardware monitor displays the pumps and valves in the live status, as well as the status of the fluid bottles.

The waste (W), air (A), and fill pump symbols (F) are illuminated in green when active. Active separator valves (Valves 1-3) are indicated in green, defective rotary valves in red. Valve status is indicated by o for open and c for closed. The fluid bottles in the hardware monitor as well as on the instrument will be illuminated in red when empty; the waste bottle will be illuminated in red when full.

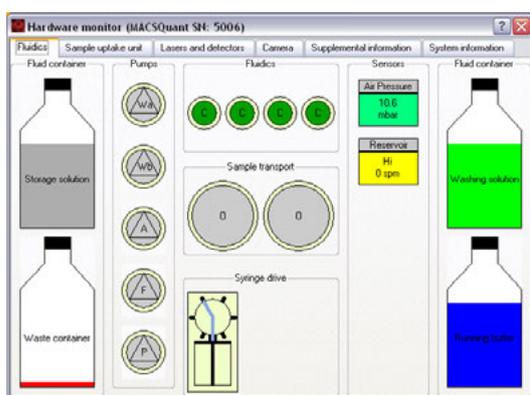
Figure 5.1: Real-time Hardware monitor of the fluidic components.

5.1.2 Sample uptake unit

The sample uptake tab indicates the status of the robotic needle arm as well as whether the MACSQuant Orbital Shaker is connected or not. The relative position of the robotic needle arm is listed in the lower left box as well as whether the Single tube rack is connected or not.

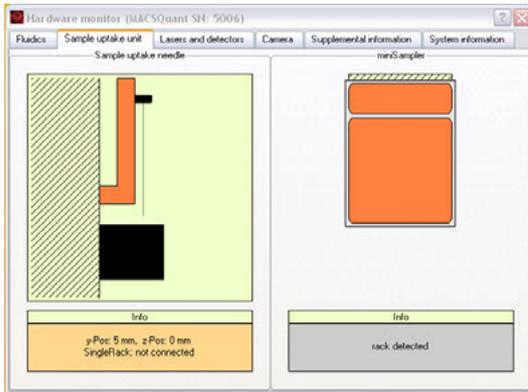
Figure 5.2: Real-time hardware monitor of the sample uptake arm.

5.1.3 Lasers and detectors

The lasers and detectors tab displays the status of the optical bench. It is possible to monitor the status of each laser fluorescence channel. The temperature, fan speed, PMT voltage, and annotated path of each laser is shown. A status overview of the optical bench is schematically represented. The bench temperature is kept between 33 °C and 37 °C. Therefore the fan speed is regulated automatically depending on the ambient room temperature and internal temperature of the MACSQuant® Instrument. Please note that the temperature for the lasers should be between 10 °C and 45 °C. In case of errors, please contact Miltenyi Biotec Technical Support.

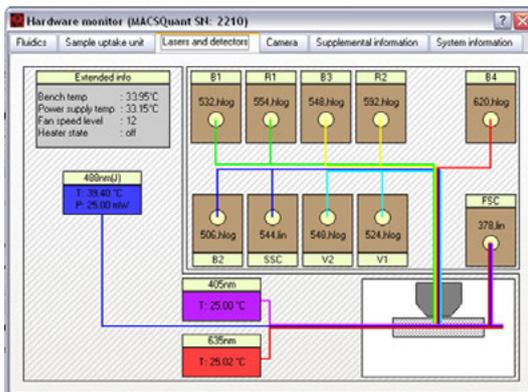
Figure 5.3: Real-time Hardware monitor of the optical bench.

6

Live Support

MACSQuant® Live Support is a real-time diagnostic service provided by Miltenyi Biotec Technical Support. Highly trained MACSQuant Specialists can be reached in real-time to assist with any queries you may have. A webcam is provided for communication with the MACSQuant Live Support, which is attached via USB connection to the MACSQuant Instrument. The webcam is installed to a free USB port at the back of the instrument and detected automatically by the software. During normal operation, the webcam should not be installed on the instrument. Please note that the MACSQuant Instrument must have network access to the internet in order to use the MACSQuant Live Support.

- 1 Select **MACSQuant Live Support** from the **Tools tab** in the **Sidebar** of the MACSQuantify Software to access remote assistance.
- 2 Complete all fields and detail any queries you may have using the **Message/Question box**.
- 3 Click **Submit** to commence MACSQuant Live Support.

7

Maintenance

7.1 Decontaminating the instrument

If the instrument has been used to process biohazardous samples, the fluidics system is to be decontaminated by running the standard shutdown sequence. It is recommended to use the appropriate disinfectant for the potential pathogen. Refer to section **Excessive debris is present in acquisition**, page 83 if you would like to perform the **Clean program** with the MACS Bleach Solution. Use tissues or swabs to decontaminate surfaces. The uptake needle of the robotic arm and the surface of the instrument shall be decontaminated upon contact with biohazardous samples.

7.2 Exchanging fluid bottles

- 1 Install one bottle at a time. Remove the bottle to be exchanged from the bottle holder
- 2 Place a new, sealed bottle into the orange bottle holder. Please note the corresponding color-coding.
- 3 Unscrew the lid of the new fluid bottle.
- 4 Remove the bottle closure from the old bottle and install it on the new bottle. Do not disconnect the color-coded tubing and sensor cable from the bottle closure. Take care not to invert the bottle closures, as this could cause liquid to enter the hydrophobic air filter.

7.3 Exchanging hydrophobic air filters

Hydrophobic air filters are attached to the bottle closures to vent the liquid bottles and the fluidic system. To avoid clogging of the filters and to prevent contamination of the liquids, exchange the hydrophobic air filters if they come into contact with liquid. To avoid clogging due to deposition of dust, exchange filters once a year. Please use only the hydrophobic air filters from Miltenyi Biotec.

- 1 To exchange the air filter of the fluidics system, remove the Single tube Rack if installed.
- 2 Open the right front cover.
- 3 Exchange the filter.
- 4 Re-mount Single tube Rack or if desired.

7.4 Cleaning the washing station

- 1 Remove the Single tube rack if needed.
- 2 Open the cover of the washing station to the left side. **Warning!** Potentially contaminated liquid may spill out of the orifice of the washing station and the tubing.
- 3 Clean the washing station by wiping it with tissue and an appropriate disinfectant, e.g. 70% ethanol, isopropyl alcohol, or 10% bleach. Finally, rinse with distilled water.

7.5 Cleaning the uptake needle

Clean the uptake needle of the robotic needle arm regularly to prevent contamination or clogging. Clean the uptake needle by wiping it with tissue and an appropriate disinfectant, e.g. 70% ethanol, isopropyl alcohol, or 10% bleach. Finally, rinse with distilled water. Move the robotic needle arm up and down to get access to the entire surface. Calibrate the robotic needle arm afterwards (refer to section **Calibrating the robotic needle arm**, page 59).

7.6 Exchange the sample needle

Before exchanging the sample needle, make sure that the needle is in its home position (that is the needle is hidden within the needle arm).



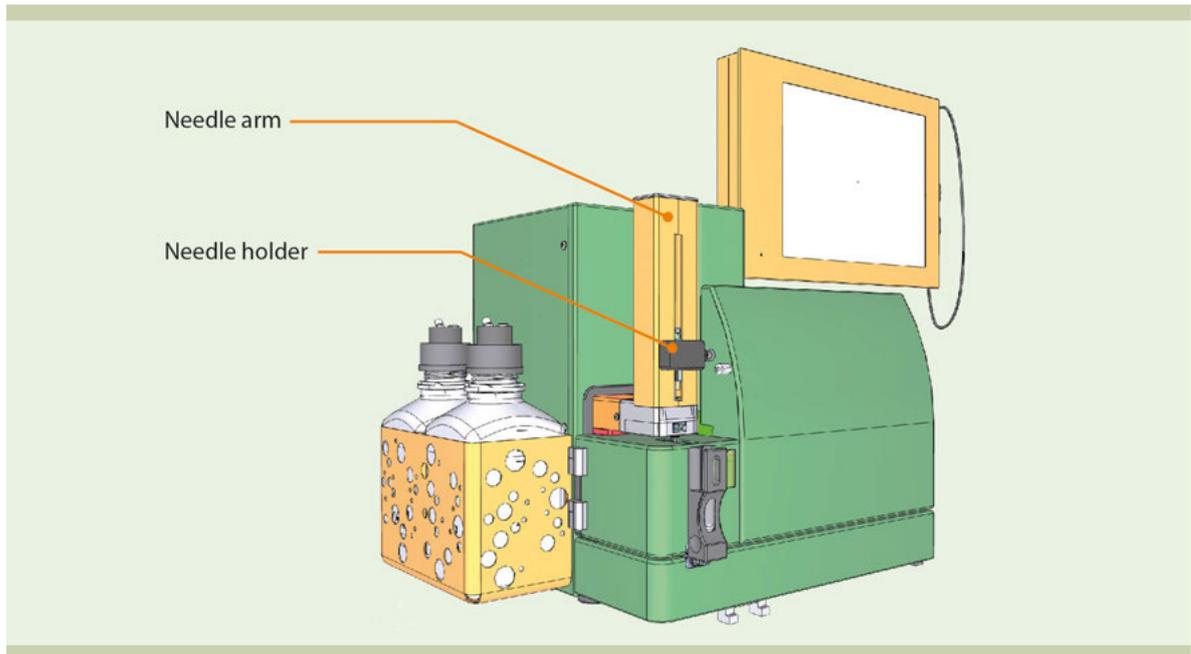
Table 7.1: Needle arm and needle holder of the MACSQuant X.


- 1 Switch off the instrument.
- 2 Remove the needle holder from the needle arm. Make sure not to bend the needle too far to the front.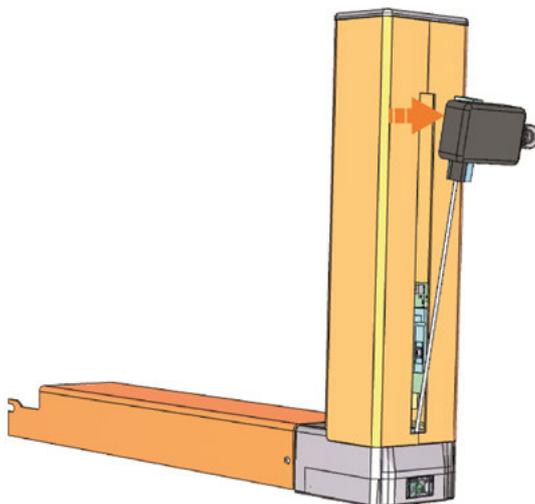


- 3 Lift the needle upwards to remove from the needle sealing.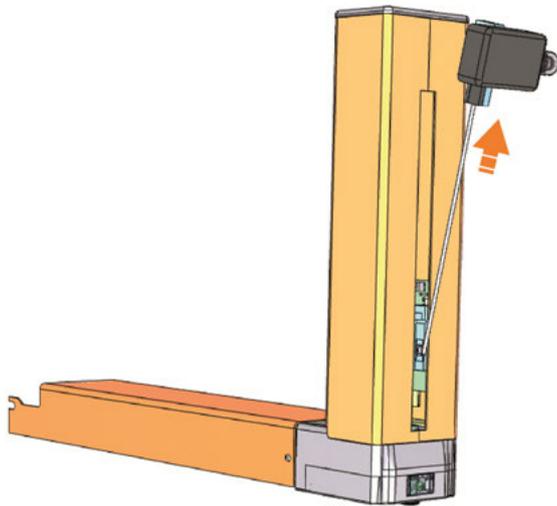


- 4 Unscrew the knurled screw of the sample tubing. Make sure not to loosen the inlay.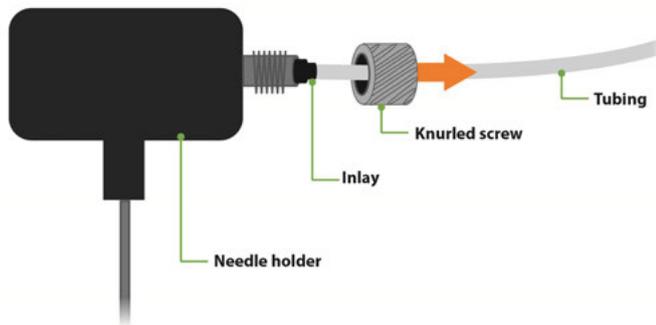

- 5 Remove the sample tubing and replace it with new tubing.
- 6 Insert the inlay into the needle holder.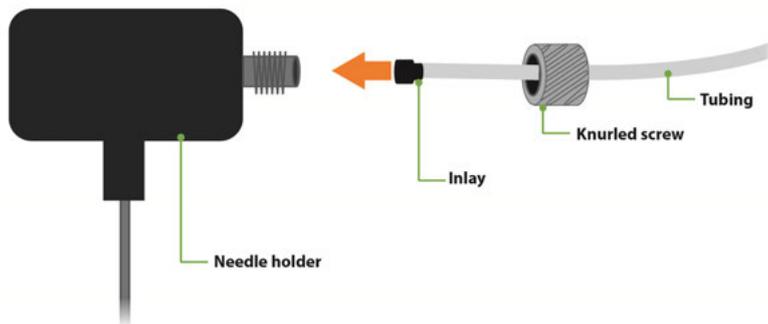

- 7 Tighten the knurled screw.

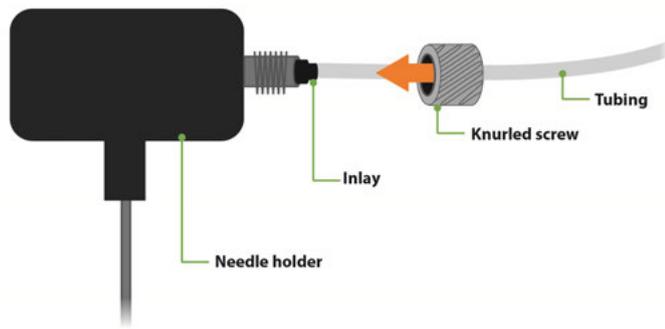- 8 Insert the needle back into the slide in the z-axis. Do not bend the needle.
- 9 Snap the needle holder back into the needle arm.
- 10 As the needle has been removed from the needle arm, its position might have been changed. Check the position and recalibrate if necessary. Mount the appropriate rack and go to **Calibrate uptake unit** to calibrate (refer to chapter **Calibration**, page 59).

7.7 Exchange the sample tubing

Exchange the sample tubing every 6 month.



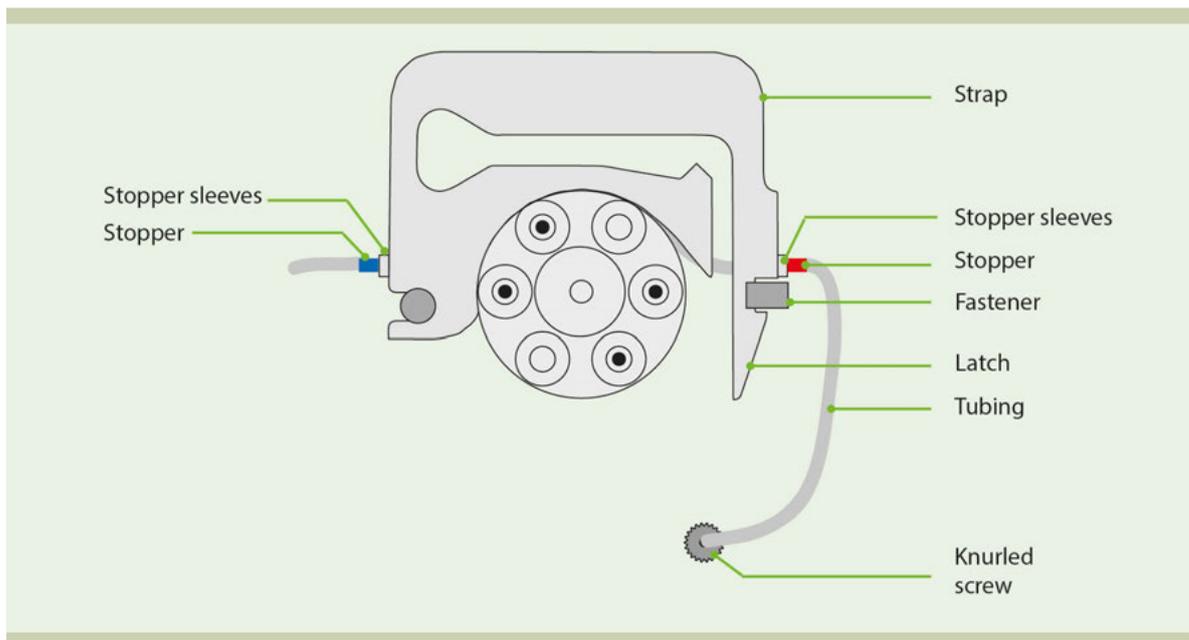
Figure 7.1: Sample tubing, mounted on tubing strap.


- 1 Switch off the instrument. Ensure that the needle is retracted completely (in the needle arm).
- 2 Open the front door.
- 3 Press the strap latch (orange arrow) to release the strap.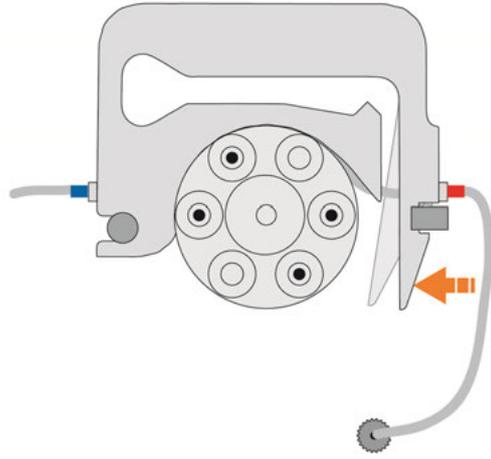


4 Lift the strap.

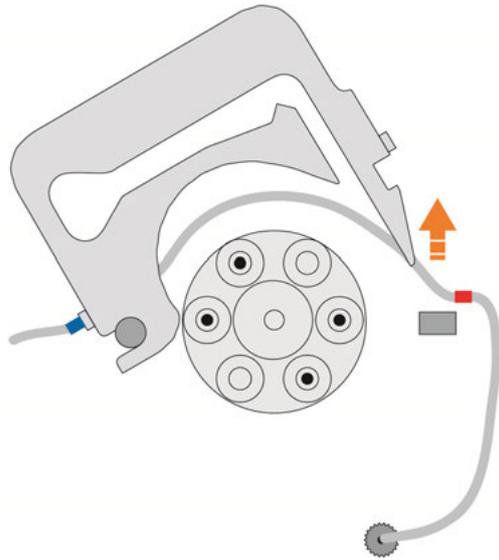
- 5 Remove the tubing strap.

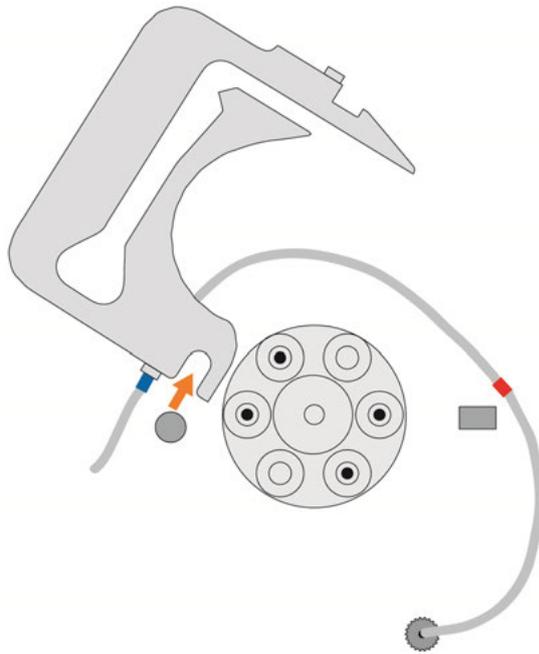- 6 Disconnect the needle from needle arm (refer to section **Exchange the sample needle**, page 74). Take care not to bend the needle.
- 7 Unscrew the knurled screw of the sample tubing. Make sure not to loosen the inlay or the stopper sleeves. Remove the sample tubing and discard.
- 8 Prepare the new sample tube: Move the knurled screw up to reveal the inlay.
- 9 Insert the tube into the needle holder.
- 10 Insert the inlay first before tightening the knurled screw.
- 11 Insert the needle back into the slide. Make sure not to bend the needle.
- 12 Connect the sample tube to the peristaltic pump unit.
- 13 Place the left side of the new tubing into the strap. Make sure that the stopper (blue) fits into the stopper sleeve.
- 14 Mount the left side of the strap, but keep it open at the right side.
- 15 Place the right side of the tubing into the strap. Make sure that the stopper (red) fits into the stopper sleeve.

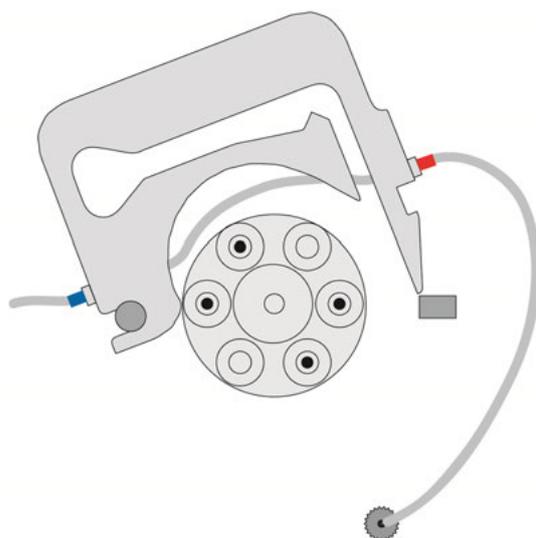16 Press the right upper corner until the strap snaps back into position.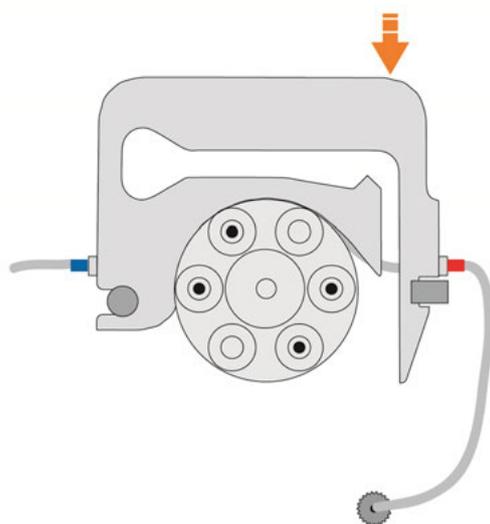

17 After exchange of the sample tubing, the system must be taught to this specific tubing in order to correctly set the tube-volume and flow rate. This process is automated. Run the ExpressMode **Volume Calibration** within the Setup Tab. The EM will guide you through this process and will take ~ 48 min. Repeat once a month.

18 The sample tubing is subject to wear, which could affect the tube-volume. In order to renew the teaching process, the ExpressMode **Tubing Volume** should be run once a week and will take ~12min.

7.7 Exchange the needle sealing

The needle sealing forms the central part of the washing station. Running buffer flows through the needle sealing perpendicular to the needle. It enables the washing of the outer wall of the needle and the buffer supply for needle and sample tubing.

The needle must be exchanged every 6 month.


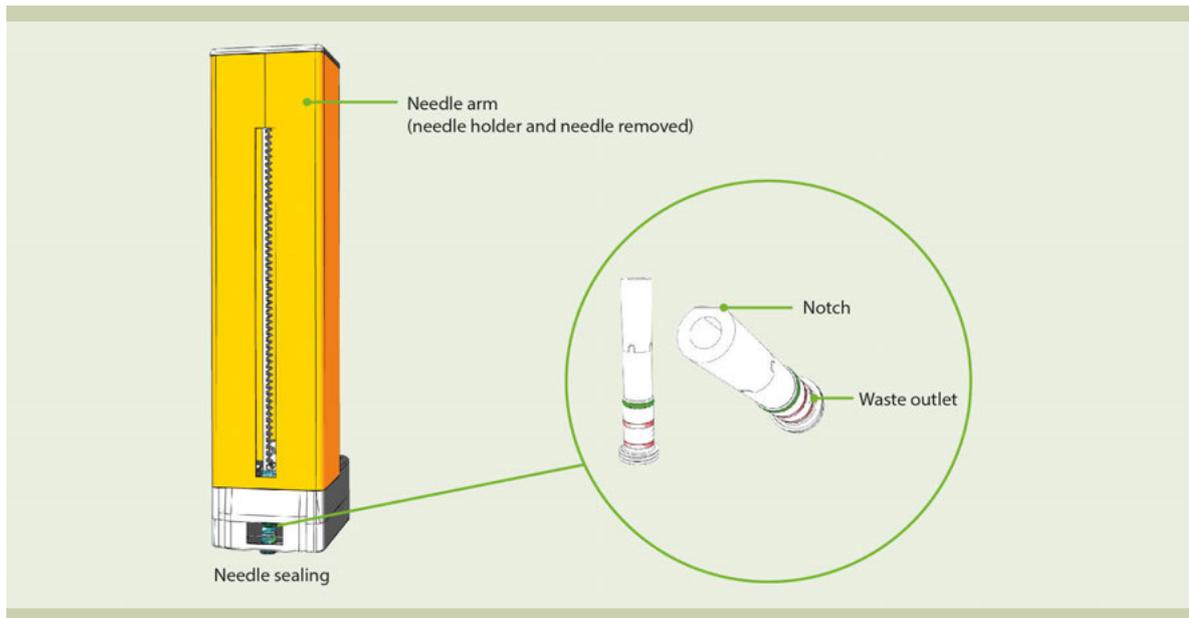
Figure 7.2: Position of the needle sealing.

- 1 Shut down the instrument.
- 2 Move needle arm to the front.
- 3 Remove needle from the needle arm as described in section **Exchange the sample needle**, page 74. You do not have to remove the sample tubing from the needle holder.
- 4 Place the sealing removal tool on the lower part of the sealing.

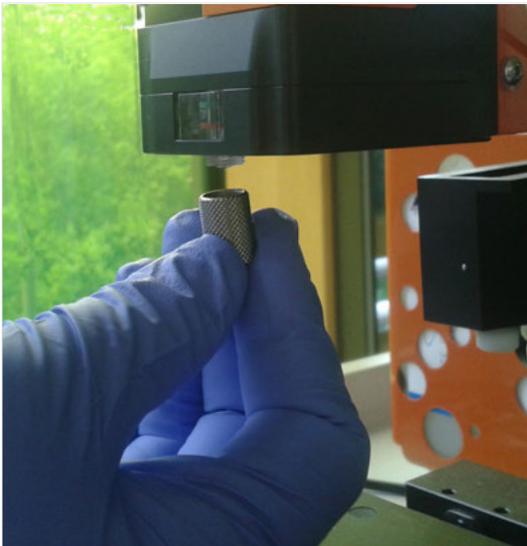- 5 Screw the removal tool counter-clockwise.
- 6 The needle sealing will move downwards until it can be pulled out.
- 7 Continue screwing to separate the sealing removal tool from the old needle sealing.
- 8 The orientation of the needle sealing is important. The top part of the sealing has a notch to indicate the correct orientation. This notch must face to the back. When viewed from the front, the top buffer-supply inlet port is facing to the left, while the lower waste outlet port is facing to the right.
- 9 Carefully push the new needle sealing from the bottom. Make sure not to tilt the needle sealing to prevent a damage of the outside sealing (colored rings).
- 10 Upon resistance the orientation of the needle sealing might have to be adjusted.
- 11 Push the needle sealing all the way in, until it snaps into final position.

- 12 Insert the needle as described in section **Exchange the sample needle**, page 74.
- 13 As the needle has been removed from the needle arm, its absolute position might have been changed. Check the position relative to the wells using the Tool **Calibrate uptake unit** (see chapter **Calibration**, page 59). Mount the appropriate rack and press **Start calibration**. If the position is still in place, confirm with **Save calibration**. Otherwise adjust the needle arm prior confirmation.

8

Troubleshooting

8.1 Syringe pump leakage

Depending on the level of use, the syringe pumps should be cleaned every 1 to 3 months. Correct overnight and long-term storage assures that no salt deposits accumulate in the syringe pumps, which could cause wear of the pump seal. The syringe pump should not run dry at any time. This damages the pump seal. In case of leakage of the syringe pump, turn off the instrument and check the connection between the syringe pump and the valve. If it is loose, tighten it. If the leakage persists, contact Miltenyi Biotec Technical Support to replace either the pump syringe or the pump seal.

8.2 Air bubbles during measurement / no event acquisition

Check for leakage within the fluidics and tubing system and perform a rinse program. Also, ensure that the sheath particle filter is devoid of air. Ensure that the needle calibration was performed correctly. Furthermore, check that the tubes connected to the bottle closures are tightly sealed. If the problem persists, contact Miltenyi Biotec Technical Support.

8.3 Excessive debris is present in acquisition

- 1 Check all fluidics to ensure that all tubes are connected properly. Ensure that the sheath particle filter is devoid of air.
- 2 Rinse or clean the system. Click the **Rinse button** in the Instrument status bar. Perform this twice.
- 3 If this does not work, try to clean the system by using the **Clean** and/or **Flush program**. To access the **Flush** program, right-click on the **Rinse button**. Click the **Clean button** to perform a system clean with MACSQuant Bleach Solution. Follow the instructions in the dialog box that appears and place the required bleach solution in the Single tube rack. The Single tube rack must be attached and selected in the **Experiment tab** in order perform a **Clean program**. The **Flush program** will perform a system backflush. This process will take approximately 16 minutes.

8.4 Touchscreen remains dark

If the touchscreen remains dark, check if the power cord is plugged in and the power supply is switched on. If the instrument is switched on. If the instrument is switched on and the touchscreen remains dark, switch off the instrument, wait five seconds and switch on again. If the instrument still does not initialize, contact Miltenyi Biotec Technical Support.

9

Technical data and specifications

The MACSQuant[®] X is labeled as a protection class I device and must be plugged into a grounded power outlet.

Please note that for USA and Canada, the main power supply cord and plug of the instrument shall comply with the following specifications: UL listed and KAM cord, minimum type SJ, minimum 18 AWG, 3 conductors. Rated for a minimum temperature of 60°C. Provided with grounding-type (NEMA 5-15P) attachment plug, rated 125 VAC, 10 A. Opposite end terminates in IEC 320-style connector, rated 125 VAC, 10A.

Supply voltage fluctuations up to +/- 10% of the nominal voltage. Transient over-voltages present on the mains supply: category II. The instrument is suitable for rated pollution degree 2.

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna,
- Increase the distance between the equipment and the receiver,
- Connect the equipment to an outlet on a circuit different from that to which the receiver is connected,
- Consult the dealer or an experienced radio/TV technician for help.

Changes or modifications of the instrument, unless expressly approved by Miltenyi Biotec, may void your authority to operate the instrument pursuant to FCC 47 CFR.

For other safety considerations, refer to the product label, or visit www.miltenyibiotec.com.

Design and specifications are subject to change without notice.

Technical data	
Model	MACSQuant X (# 130-105-100)
Color	Green/orange
Footprint (w × d)	669 mm × 368 mm (26.3×14.5")
Footprint with MACSQuant X Orbital Shaker (w × d)	814 mm × 473 mm (32×18.6")
Height (with MACSQuant X Orbital Shaker)	392.5-582 mm (adjustable touchscreen)
Weight	50 kg (110 lbs)
Input voltage	100-240 VAC, 50/60 Hz
Power consumption	450 W
Fuses	2 x 8 AT, 250 V
RS232 Interface (labeled "RS232/AUX") Not in use	Pin1, 4, 6-9: NC Pin 2: RXD Pin 3: TXD Pin 5: GND
AC Output (labeled "Bottle Sensor")	Pin1-5: Input Pin 6-15: GND Pin 9-13: Input/Output 5 VAC/10kΩ
RS232 Interface + DC-Output (labeled "RS232/BCR")	Pin1: Input/Output Pin 2:TXD Pin 3: RXD Pin 4,6: NC Pin 5: GND Pin 7, 8: Shorted Pin 9: 5 VDC/2.6 A
CAN Bus + DC-Output (labeled "External CAN")	Pin 1, 4, 8: NC Pin 2: CAN-L Pin 3, 6: GND Pin 5,9: 24 VDC/1.85 A Pin 7: CAN-H
CAN Bus + DC-Output (labeled "Power CAN")	Pin 1: CAN-L Pin 2: CAN-H Pin 3: GND Pin 5: 24 V DC, 1.1 A A1: 24 V DC, 5 A A2: GND
RS232 Interface (labeled "COM 1")	Pin 1: DCD Pin 2: RxD Pin 3: TxD Pin 4: DTR Pin 5: GND Pin 6: DSR Pin 7: RTS Pin 8: CTS

Technical data	
	Pin 9: RI
4 x USB 2.0 ports (labeled "USB 1-4")	Pin 1: +5 V Pin 2: - Pin 3: + Pin 4: GND
VGA port (labeled "VGA")	Video graphics array analog interface supports up to 2048 × 1536 @75 Hz Pin 1: RED Pin 2: GREEN Pin 3: BLUE Pin 4: ID2/RES Pin 5-8, 10: GND Pin 9: KEY/PWR Pin 11: ID0/RES Pin 12: ID1/SDA Pin 13: HSync Pin 14: VSync Pin 15: ID3/SCL
2 × HDMI ports	2×High-Definition Multimedia Interfaces support up to 1920×1200 @60 Hz Pin 1-3: TMDS Data2 Pin 4-6: TMDS Data1 Pin 7-9: TMDS Data0 Pin 10-12: TMDS Clock Pin 13: CEC Pin 14: Reserved Pin 15: SCL Pin 16: SDA Pin 17: GND Pin 18: +5 V/50 mA Pin 19: Hot Plug detect
2 x Ethernet ports (GbE) (labeled "LAN 1" and "LAN 2")	Pin1: TXD+ Pin 2:TXD+ Pin 3: RXD+ Pin 4,5,7,8: NC Pin 6: RXD-
Audio ports	Green: headphones Pink: microphone
CF Card Slot	The inserted card contains the operating system of the MACSQuant Instrument (Microsoft Windows 7 with MACSQuantify Software, read-only mode)
RAM	8 GB DDR4 (SO-DIMM, 1066 MHz)
HDD	2.5" automotive & extended temperature range
Monitor	15.6" HD touchscreen, height-adjustable
Working temperature	+15 °C to +25 °C
Humidity	10% to 90% relative humidity, non-condensing
Altitude	Max. 2000 m
Emission sound pressure level at workstation	< 70 dB(A)

Table 9.1: Technical specifications of the MACSQuant X.

9.1 EC Declaration of conformity

This declaration of conformity is issued under the sole responsibility of the manufacturer:

Miltenyi Biotec GmbH
Friedrich-Ebert-Str. 68
51429 Bergisch Gladbach
Germany

The declaration of conformity refers to the machinery identified as follows:

Description: Laboratory equipment

Model: MACSQuant® X

The machinery complies with all essential requirements of the following directives

- 2014/30/EU Electromagnetic compatibility
- 2006/42/EC Machinery

The machinery is in conformity with the following harmonized standards:

- EN 61010-2-081
- EN 60825-1
- EN 61326-1

9.1.1 Additional international conformity

The MACSQuant® X is in conformity with:

- UL 61010-1
- CAN/CSA-22.2 No. 61010-1
- IEC 61010-2-081

Parameter	Specification																						
Minimum sample volume	20 µL (recommended)																						
Scatter resolution	Scatter performance is optimized for resolving human peripheral blood lymphocytes, monocytes, and granulocytes																						
Sample flow rate	25-100 µl/min plus automated flow rate to maintain 500, 1000, or 2000 event/s/second																						
Maximal event rate	Up to 15,000 events/second																						
Lasers	488 nm, 404 nm, 640 nm																						
Fluorescence sensitivity (MESF)	FITC < 150; PE < 100; APC < 150																						
Absolute cell count performance	Accuracy +/- 5%; reproducibility (CV) <5%																						
Sample carryover	Typical value of 0.01%																						
Emission detectors	<table border="1"> <thead> <tr> <th>Channel</th> <th>Filter</th> </tr> </thead> <tbody> <tr> <td>FSC</td> <td>488/10 nm</td> </tr> <tr> <td>SSC</td> <td>488/10 nm</td> </tr> <tr> <td>V1</td> <td>450/50 nm</td> </tr> <tr> <td>V2</td> <td>525/50 nm</td> </tr> <tr> <td>B1</td> <td>525/50 nm</td> </tr> <tr> <td>B2</td> <td>585/40 nm</td> </tr> <tr> <td>B3</td> <td>655-730 nm</td> </tr> <tr> <td>B4</td> <td>750 nm LP</td> </tr> <tr> <td>R1</td> <td>655-730 nm</td> </tr> <tr> <td>R2</td> <td>750 nm LP</td> </tr> </tbody> </table>	Channel	Filter	FSC	488/10 nm	SSC	488/10 nm	V1	450/50 nm	V2	525/50 nm	B1	525/50 nm	B2	585/40 nm	B3	655-730 nm	B4	750 nm LP	R1	655-730 nm	R2	750 nm LP
Channel	Filter																						
FSC	488/10 nm																						
SSC	488/10 nm																						
V1	450/50 nm																						
V2	525/50 nm																						
B1	525/50 nm																						
B2	585/40 nm																						
B3	655-730 nm																						
B4	750 nm LP																						
R1	655-730 nm																						
R2	750 nm LP																						

Table 9.2: Operation details of the MACSQuant® X.

10

Technical support

For technical support, please contact your local Miltenyi Biotec representative or the Technical Support team at Miltenyi Biotec headquarters:

Miltenyi Biotec GmbH
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macstec@miltenyibiotec.de

Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

11

Limited warranty

Except as stated in a specific warranty statement, which may accompany this product, or unless otherwise agreed in writing by an authorized representative of Miltenyi Biotec, Miltenyi Biotec's warranty, if any, with respect to this product is subject to the terms and conditions of sale of the company within the Miltenyi Biotec group, which supplied the product. Terms and conditions of sale may vary by country and region. Nothing in this document should be construed as constituting an additional warranty.

Miltenyi Biotec's warranty for this product only covers product issues caused by defects in material or workmanship during ordinary use; it does not cover product issues caused by any other reason, including but not limited to product issues due to use in a manner other than specifically described in this manual, for example: inappropriate or improper use; incorrect assembly or installation by an operator or a third party; reasonable wear-and-tear; negligent or incorrect handling, servicing, or maintenance; non-adherence to the operating instructions; unauthorized modification of or to any part of this product; or use of inappropriate accessories or work materials.

Miltenyi Biotec's warranty does not cover products sold AS IS or WITH ALL FAULTS, or which had its serial number defaced, altered or removed, or any consumables, or parts identified as being supplied by a third party. Miltenyi Biotec must be informed immediately, if a claim is made under such warranty. If a material or manufacturing defect occurs within the warranty period, Miltenyi Biotec will take the appropriate steps to restore the full usability of your product.

Limitation on damages: Miltenyi Biotec shall not be liable for any incidental or consequential damages for breach of any express or implied warranty or condition on this product.

Some states or jurisdictions do not allow the exclusion or limitation of incidental or consequential damages, so the above limitations or exclusions may not apply to you. This warranty statement gives you specific legal rights and you may have other rights, which may vary from country to country or jurisdiction to jurisdiction.

Miltenyi Biotec

Germany/Austria/ Switzerland

Miltenyi Biotec GmbH
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macs@miltenyibiotec.de

USA/Canada

Miltenyi Biotec Inc.
2303 Lindbergh Street
Auburn, CA 95602, USA
Phone 800 FOR MACS
Phone +1 530 888 8871
Fax +1 877 591 1060
macs@miltenyibiotec.com

Australia

Miltenyi Biotec
Australia Pty. Ltd.
Unit 16A, 2 Eden Park Drive
Macquarie Park NSW 2113
Australia
Phone +61 2 8877 7400
Fax +61 2 9889 5044
macs@miltenyibiotec.com.au

www.miltenyibiotec.com

Benelux

Miltenyi Biotec B.V.
Schipholweg 68 H
2316 XE Leiden
The Netherlands
macs@miltenyibiotec.nl

Customer service

The Netherlands
Phone 0800 4020120
Fax 0800 4020100

Customer service Belgium

Phone 0800 94016
Fax 0800 99626

Customer service Luxembourg

Phone 800 24971
Fax 800 24984

China

Miltenyi Biotec Technology &
Trading (Shanghai) Co., Ltd.
Rooms 2303 and 2309
No. 319, Xianxia Road
Changning District
200051 Shanghai 200051, P.R. China
Phone +86 21 62351005
Fax +86 21 62350953
macs@miltenyibiotec.com.cn

France

Miltenyi Biotec SAS
10 rue Mercœur
75011 Paris, France
Phone +33 1 56 98 16 16
Fax +33 1 56 98 16 17
macs@miltenyibiotec.fr

Italy

Miltenyi Biotec S.r.l.
Via Paolo Nanni Costa, 30
40133 Bologna
Italy
Phone +39 051 6 460 420
Fax +39 051 6 460 499
macs@miltenyibiotec.it

Japan

Miltenyi Biotec K.K.
Nittsu-Eitai Building 5F
16-10 Fuyuki, Koto-ku,
Tokyo 135-0041, Japan
Phone +81 3 5646 8910
Fax +81 3 5646 8911
macs@miltenyibiotec.jp

Nordics and Baltics

Miltenyi Biotec Norden AB
Scheelevägen 17
223 70 Lund
Sweden
macs@miltenyibiotec.se

Customer service Sweden

Phone 0200-111 800

Fax 046-280 72 99

Customer service Denmark

Phone 80 20 30 10

Fax +46 46 280 72 99

Customer service

**Norway, Finland, Iceland
and Baltic countries**

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Miltenyi Biotec provides products and services worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for therapeutic or diagnostic use. MACS, MACSQuant and MACSQuantify are registered trademarks or trademarks of Miltenyi Biotec GmbH and/or its affiliates in various countries worldwide. All other trademarks mentioned in this publication are the property of their respective owners and are used for identification purposes only.

Copyright © 2017 Miltenyi Biotec GmbH and/or its affiliates. All rights reserved.

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Miltenyi Biotec Korea Co., Ltd
Miltenyi Biotec Ltd.